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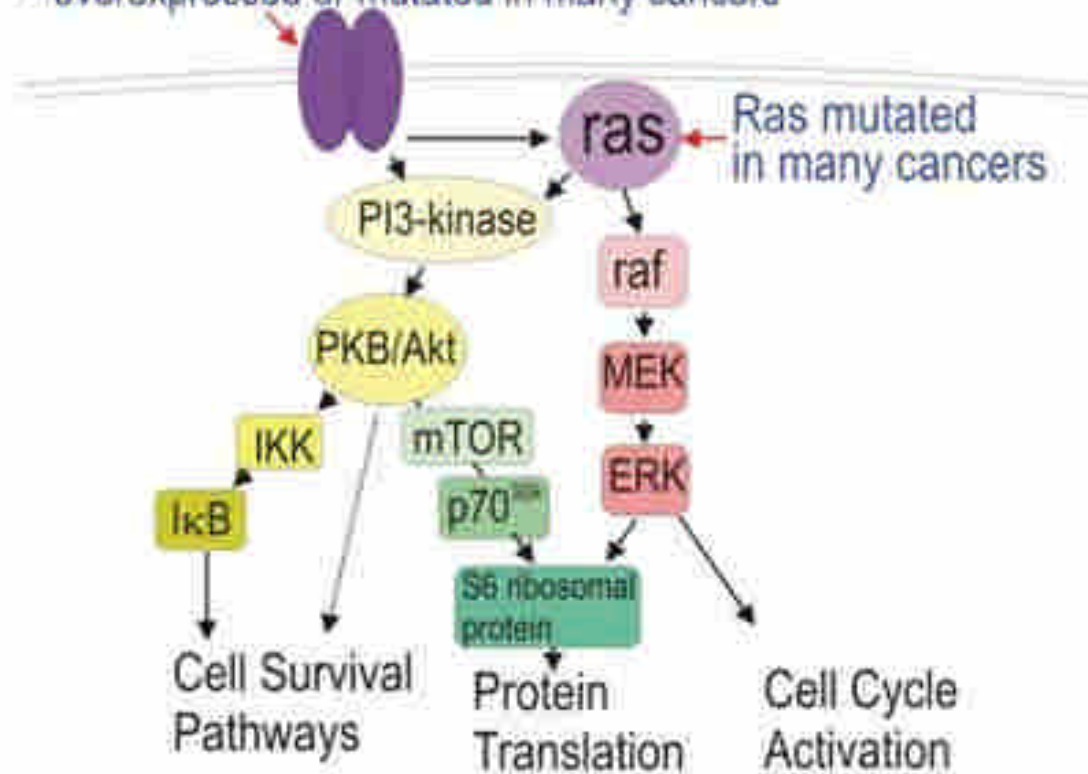
Helping all people
live healthy lives

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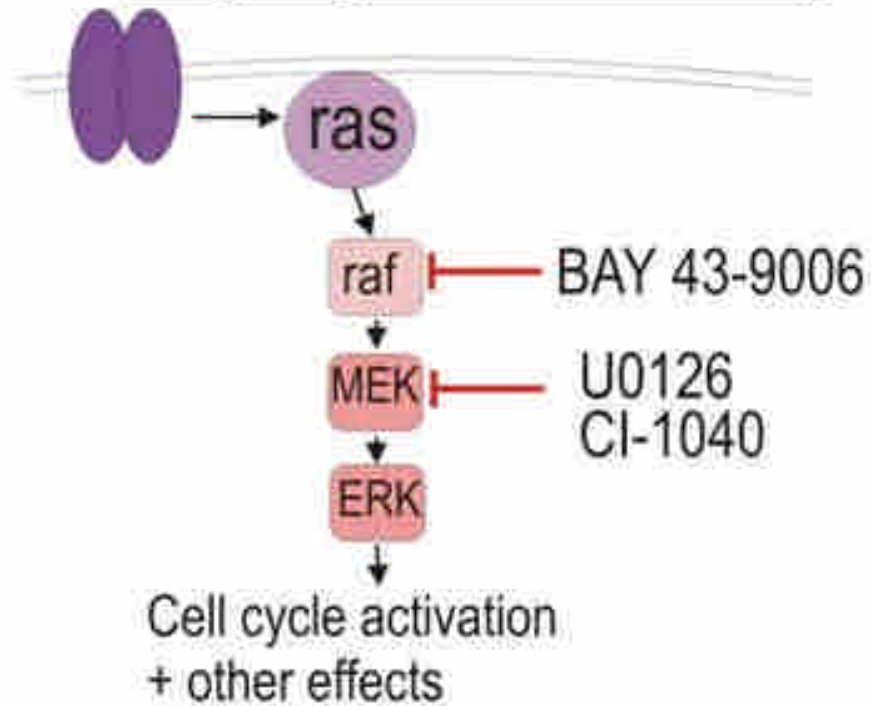
Receptor tyrosine kinases -
e.g. EGFR, Her2

- overexpressed or mutated in many cancers

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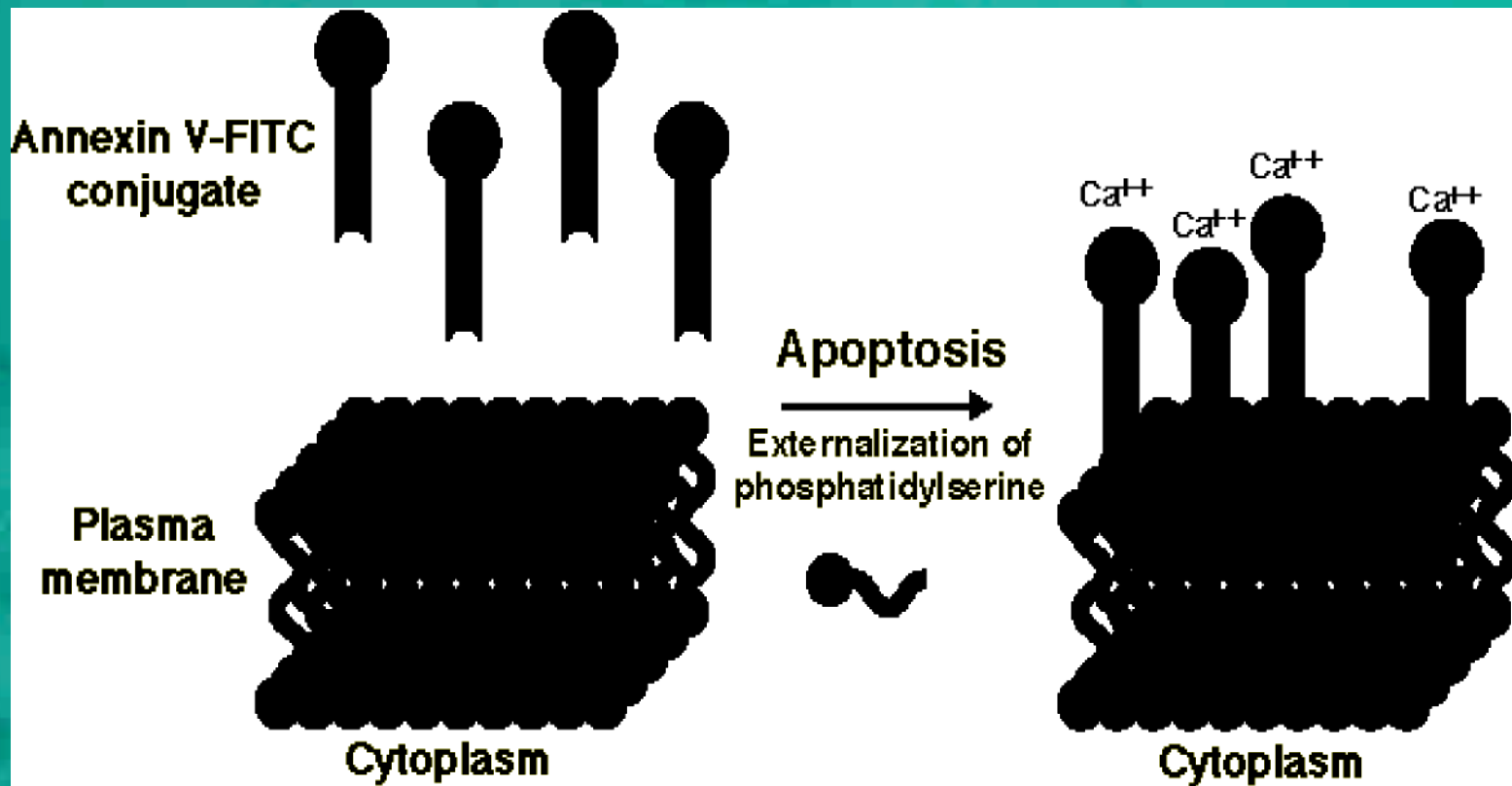
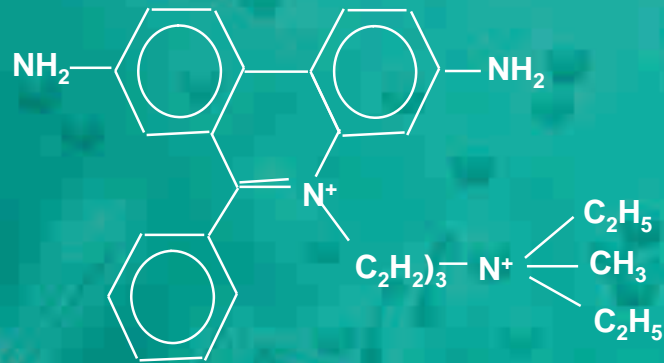
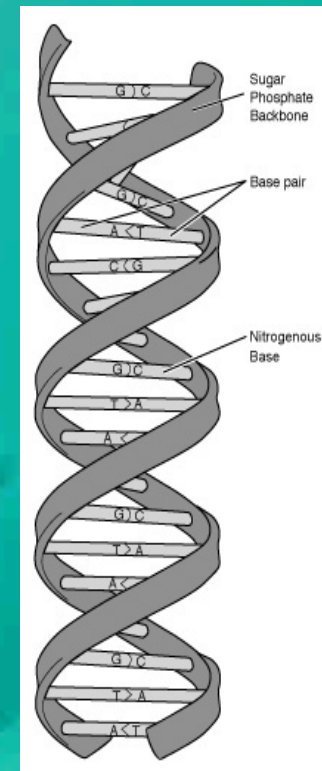
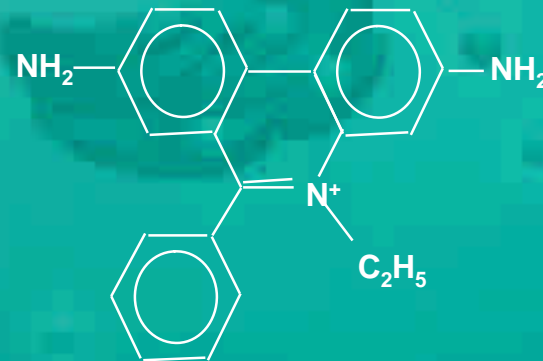
Helping all people
live healthy lives

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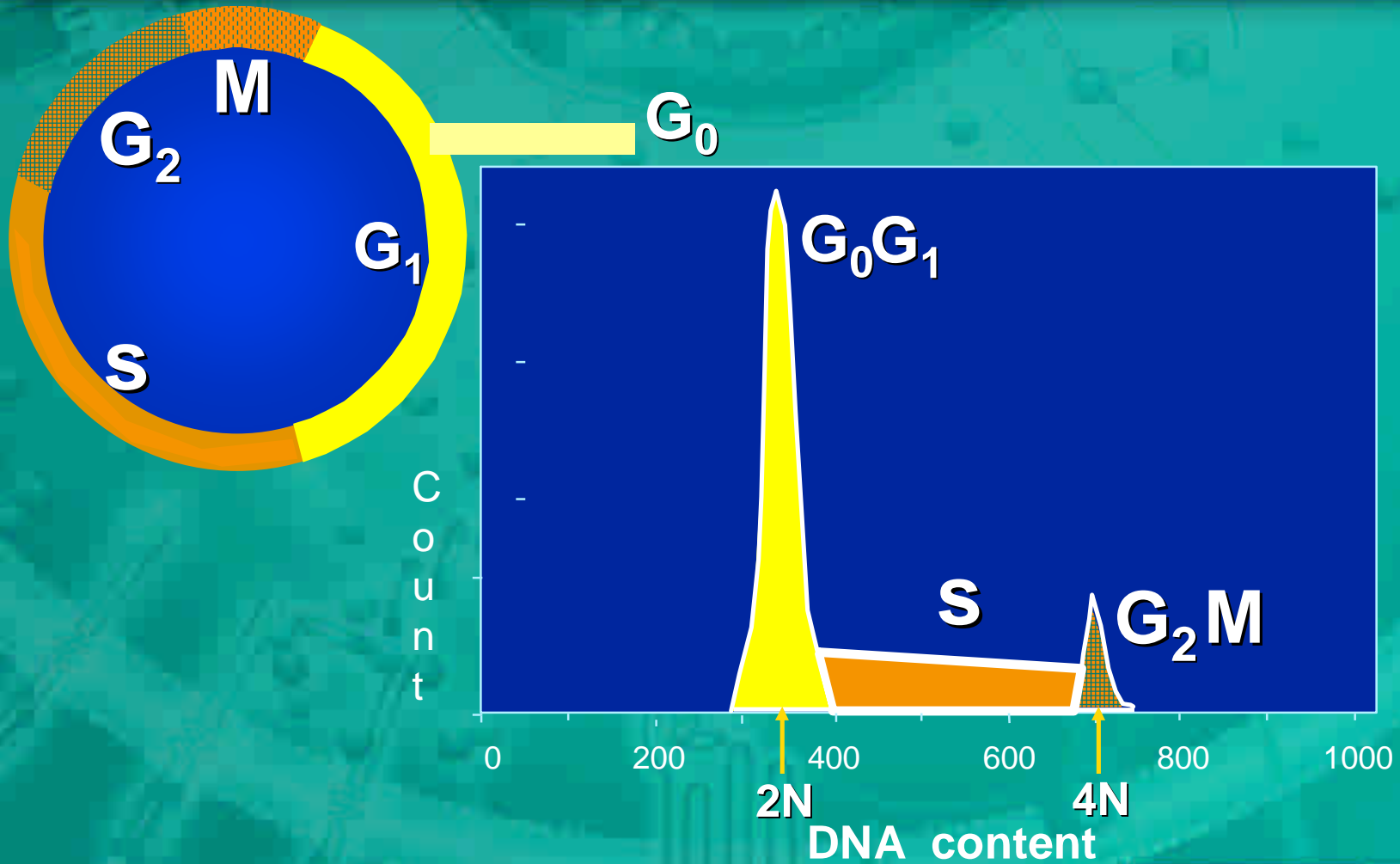
Propidium

Ethidium

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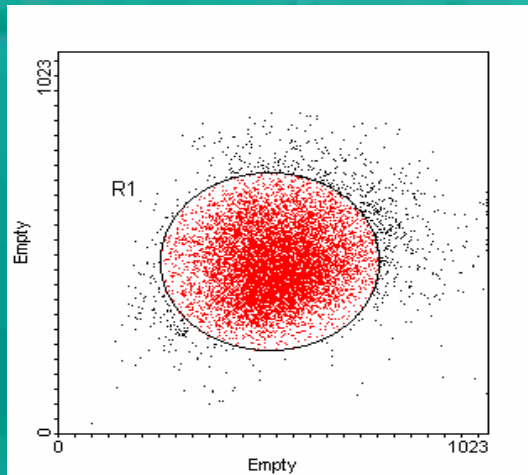
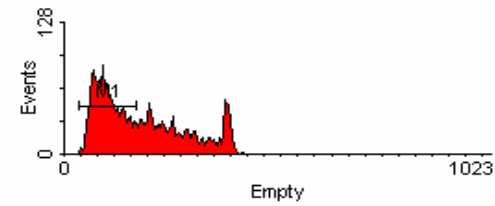
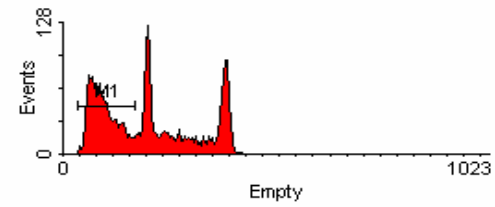
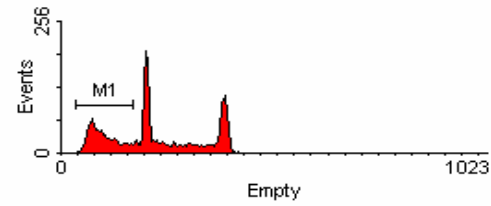
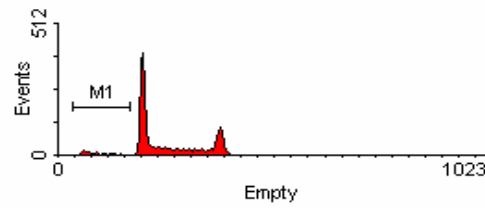
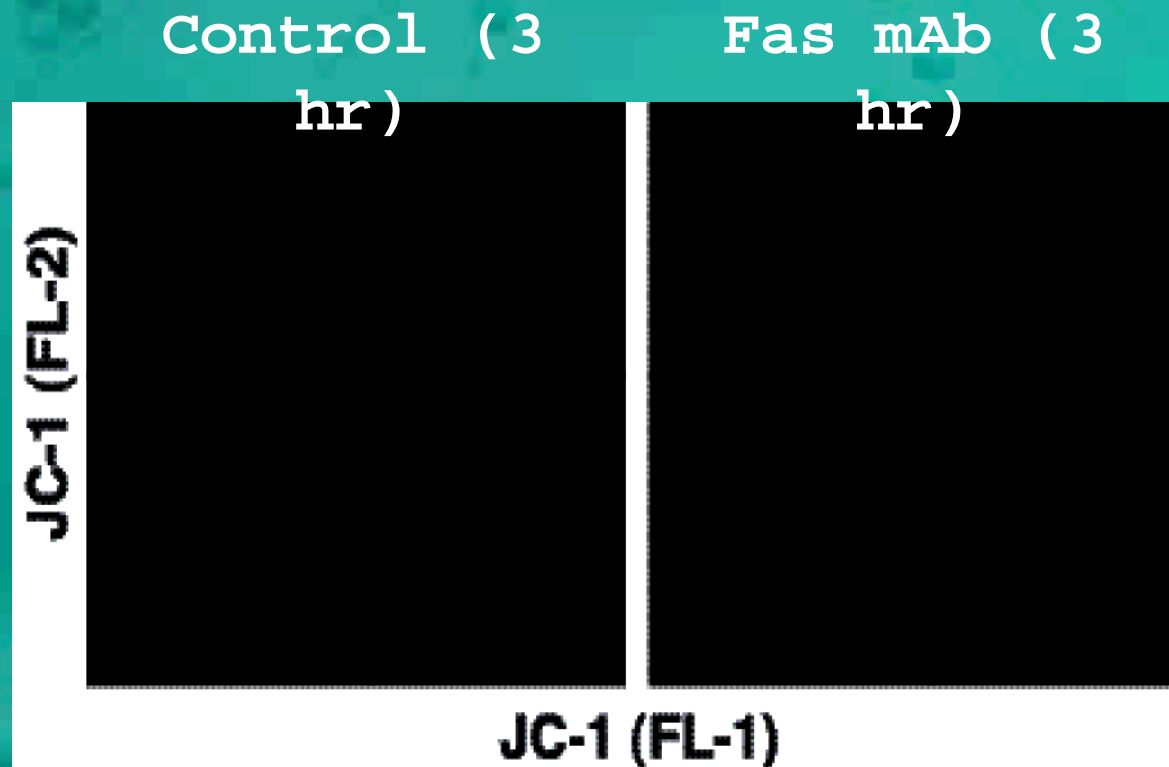
Helping all people
live healthy lives

Sub G1

Helping all people
live healthy lives

Mitochondria Membrane Potential

Helping all people
live healthy lives

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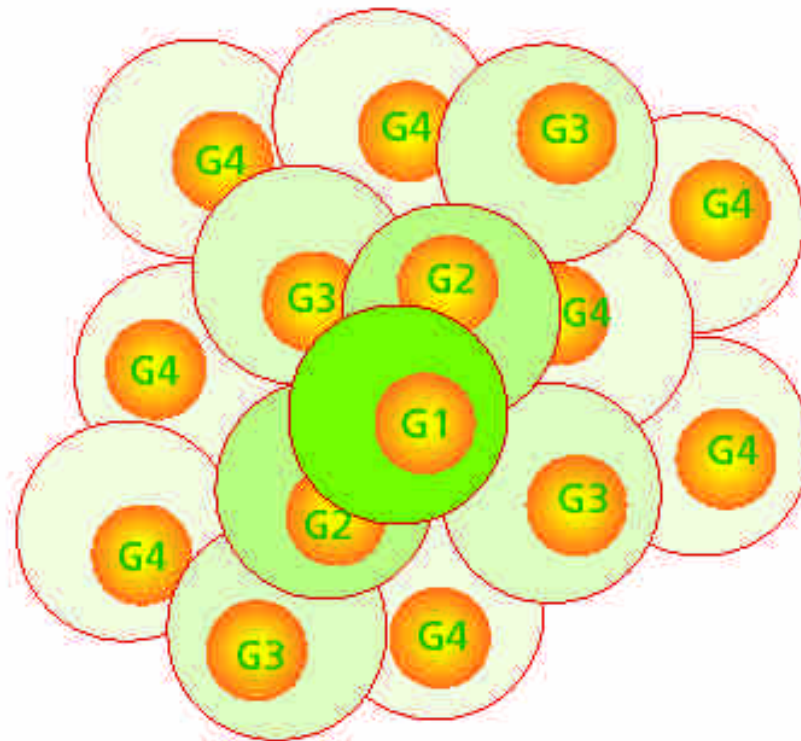
Figure 1. Upon cell division, CFSE is distributed uniformly between daughter cells.

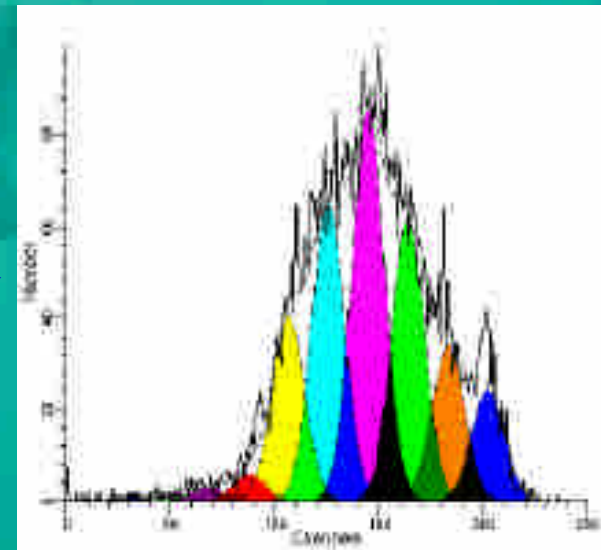
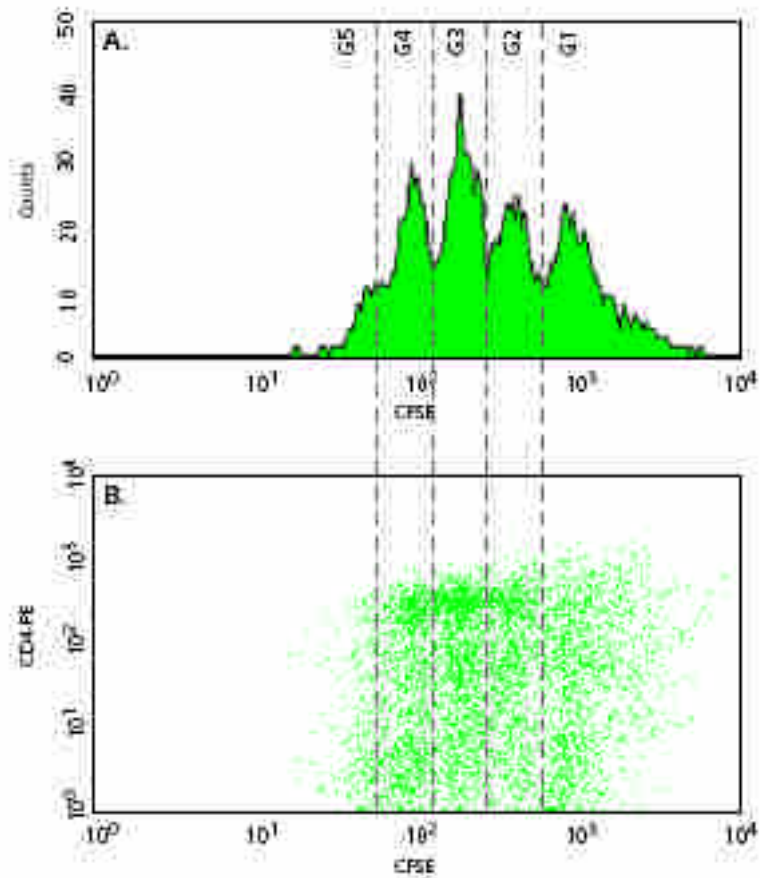
Figure 2. Profile of CFSE-labeled, PHA-stimulated (72 hrs) peripheral blood mononuclear cells (PBMCs) analyzed by flow cytometry

Helping all people live healthy lives

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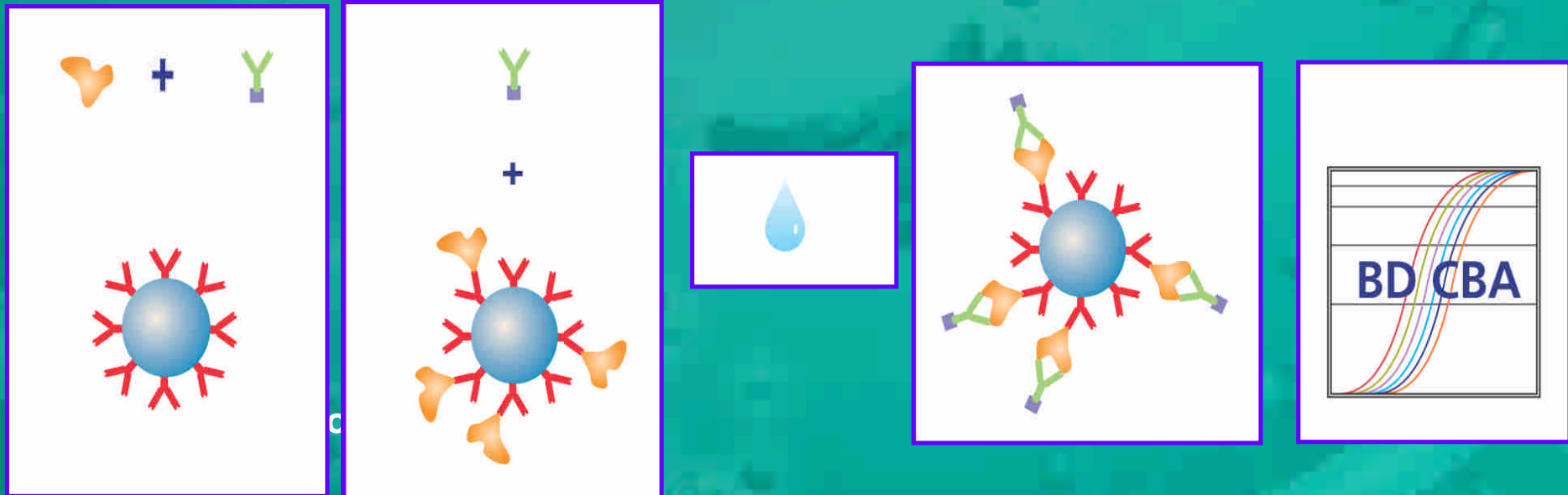
Helping all people
live healthy lives

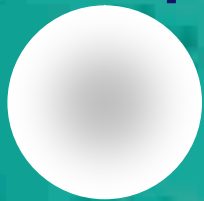
Cytometric Beads Array (CBA)

Helping all people
live healthy lives

Beads Provide a Flexible Platform

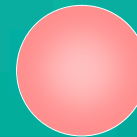
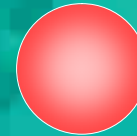
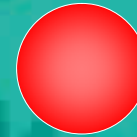
Beads provide an expandable assay platform for use with a flow cytometer

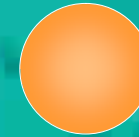
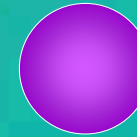
Multiple sizes

Different intensities*

Different colors with different intensities

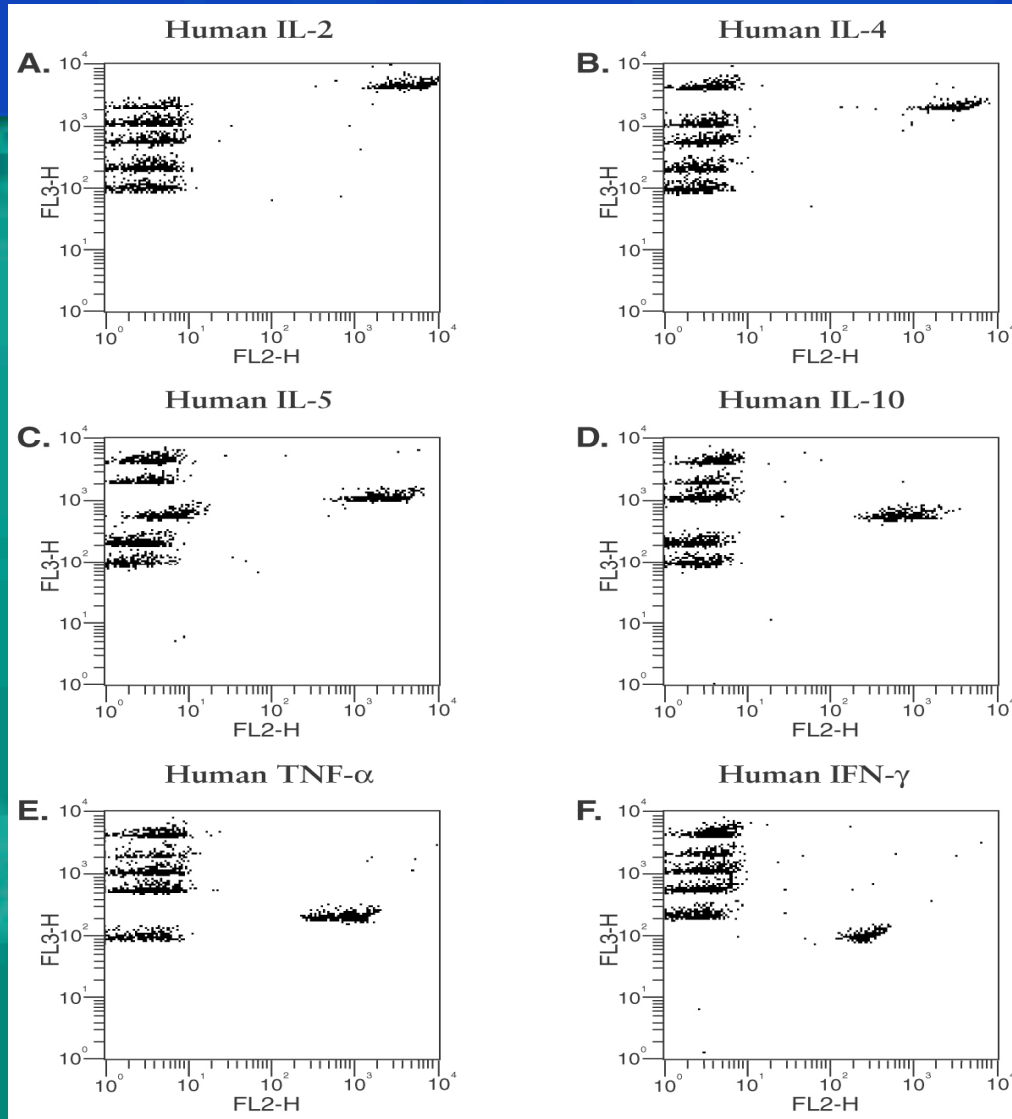
Helping all people live healthy lives

Advantages of Bead-Based Immunoassays

- Analyze multiple analytes simultaneously
- Reduced sample volume requirements
- Reduced hands-on time by parallel analysis of samples
- Wide dynamic range of fluorescence detection (requires fewer sample dilutions)

Helping all people
live healthy lives

Proteins Measured

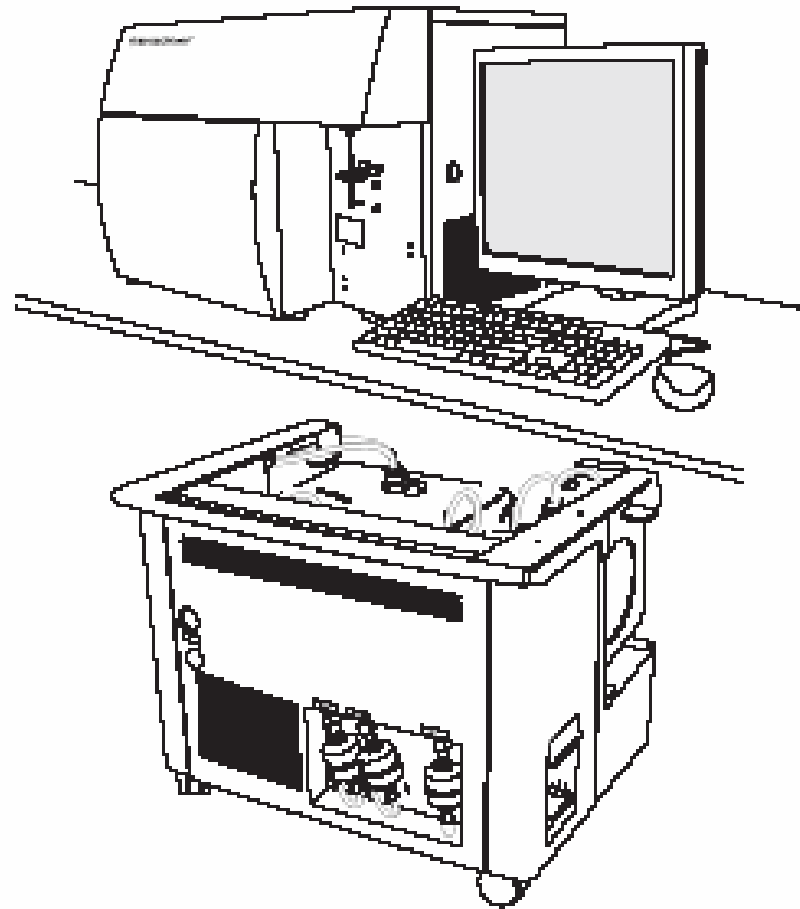
- A. Interleukin (IL)-2
- B. IL-4
- C. IL-5
- D. IL-10
- E. Tumor Necrosis Factor- α
- F. Interferon- γ

Helping all people
live healthy lives

BD FACSCanto™

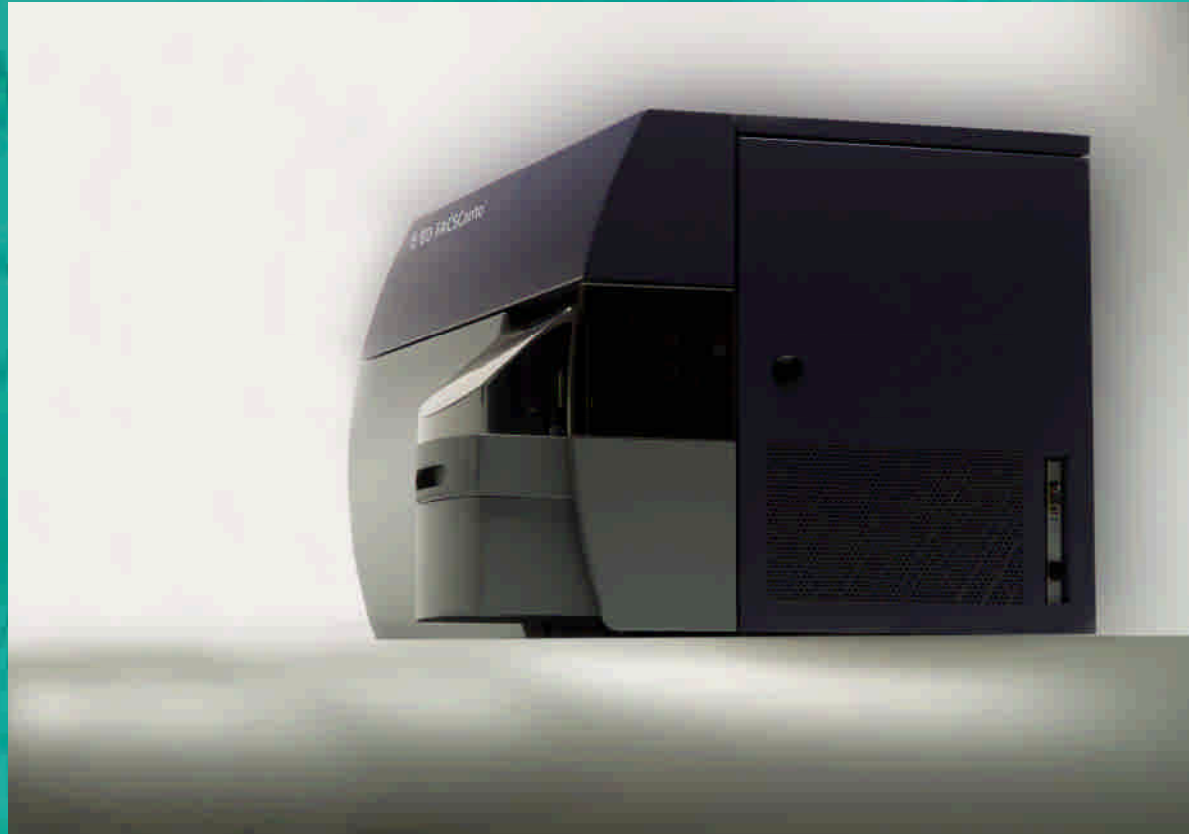
Figure 1-1 BD FACSCanto system

Helping all people
live healthy lives

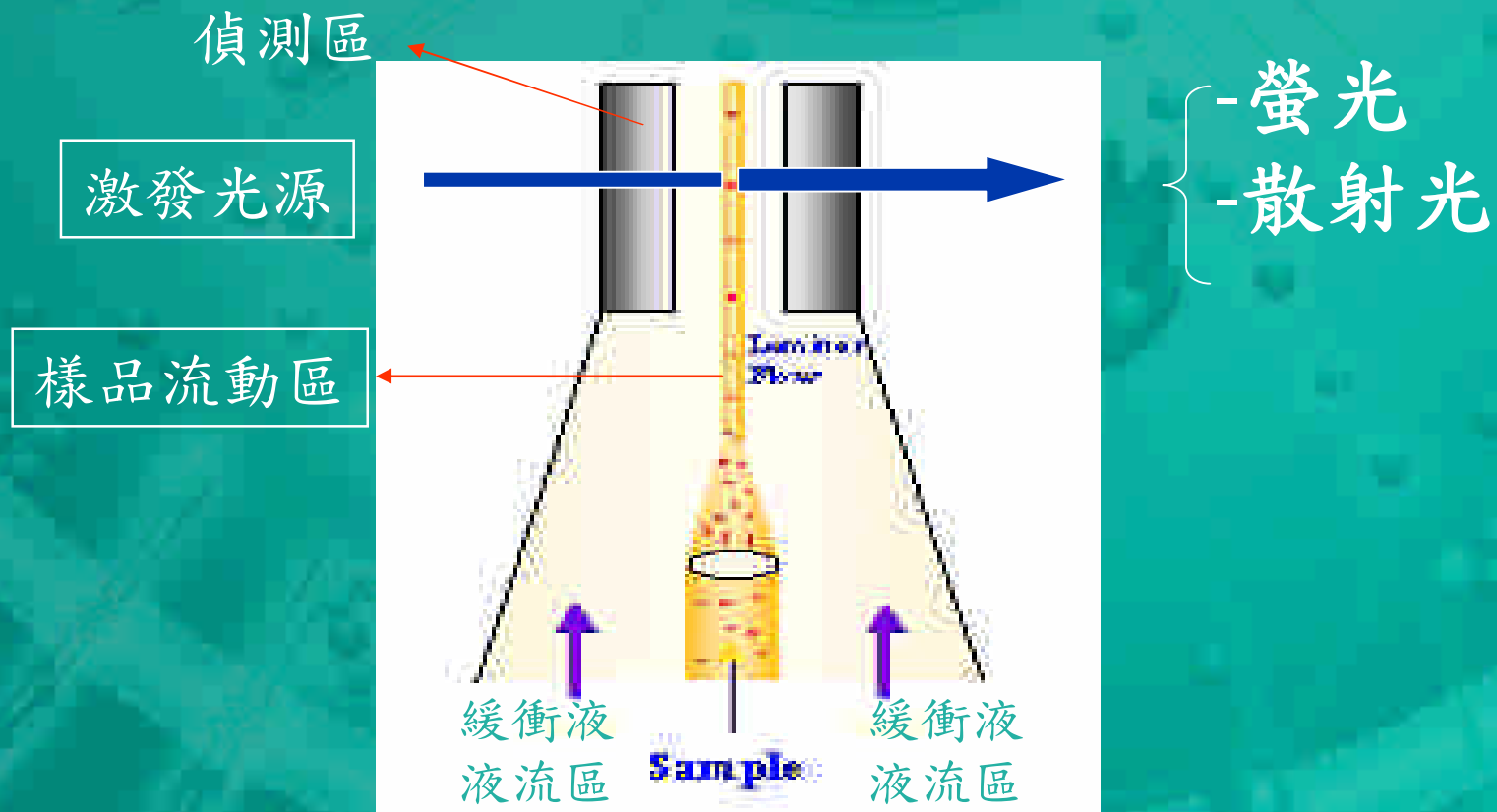
Helping all people
live healthy lives

Helping all people
live healthy lives

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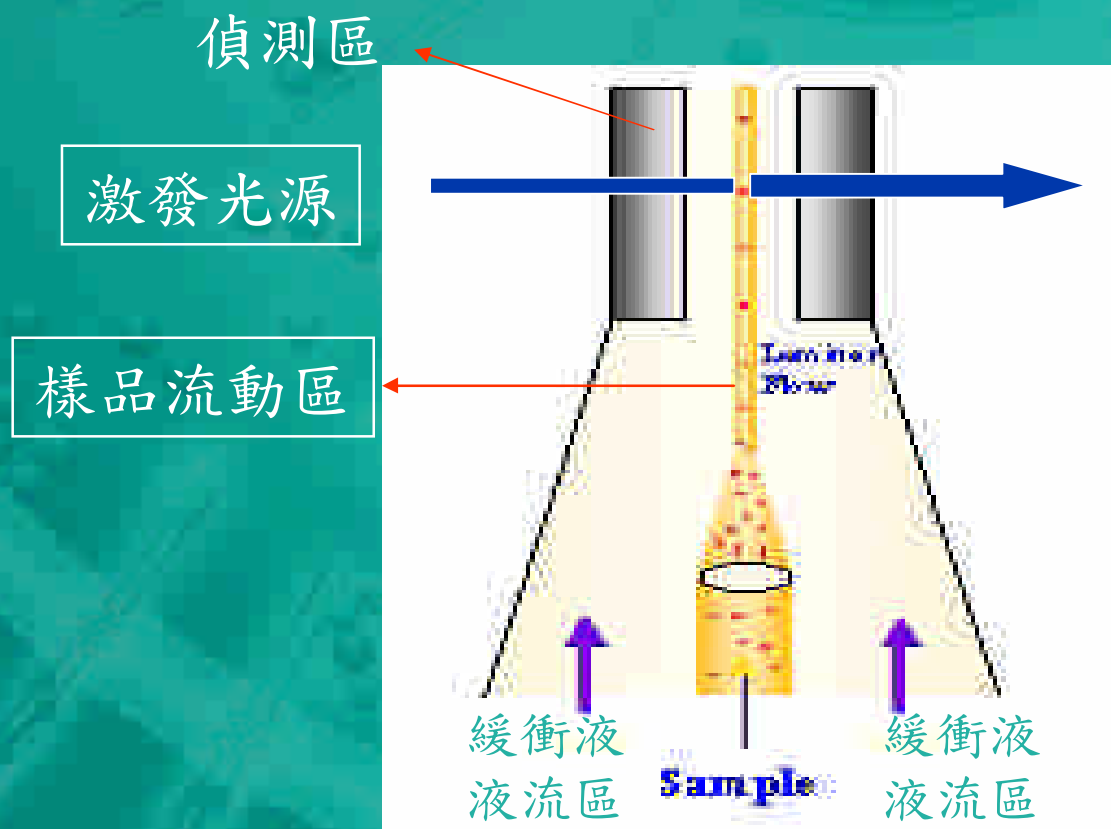
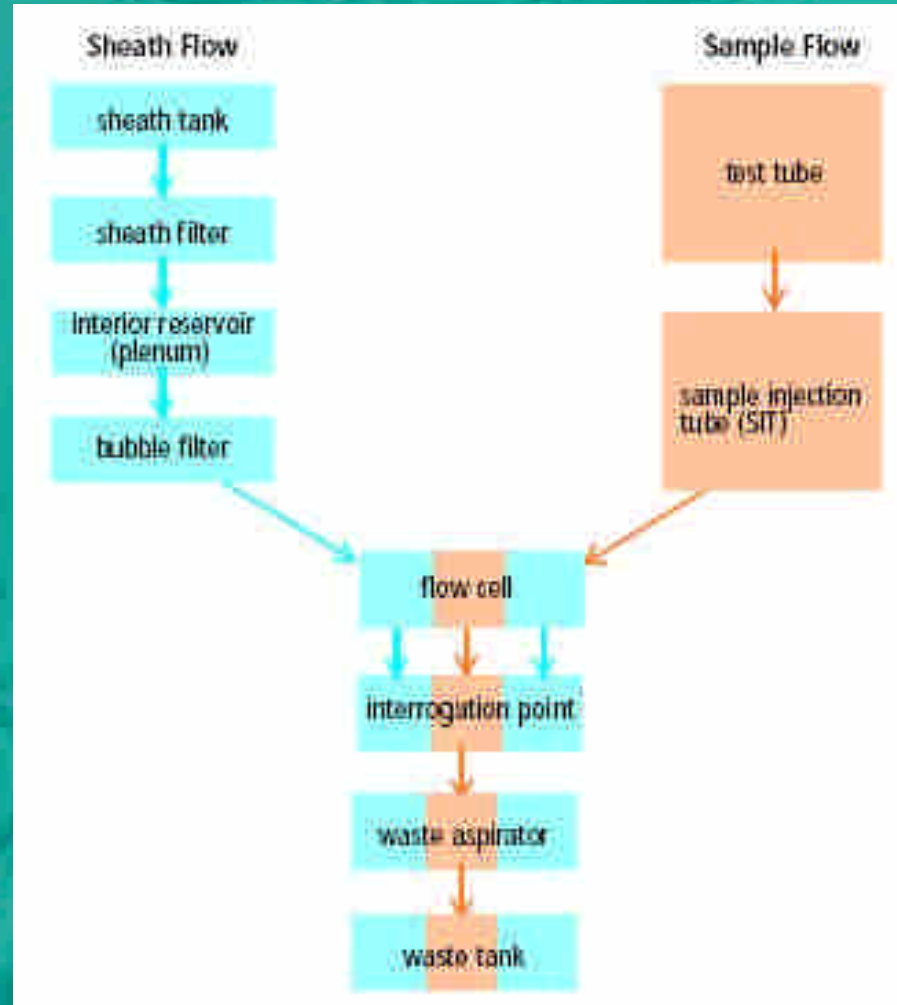
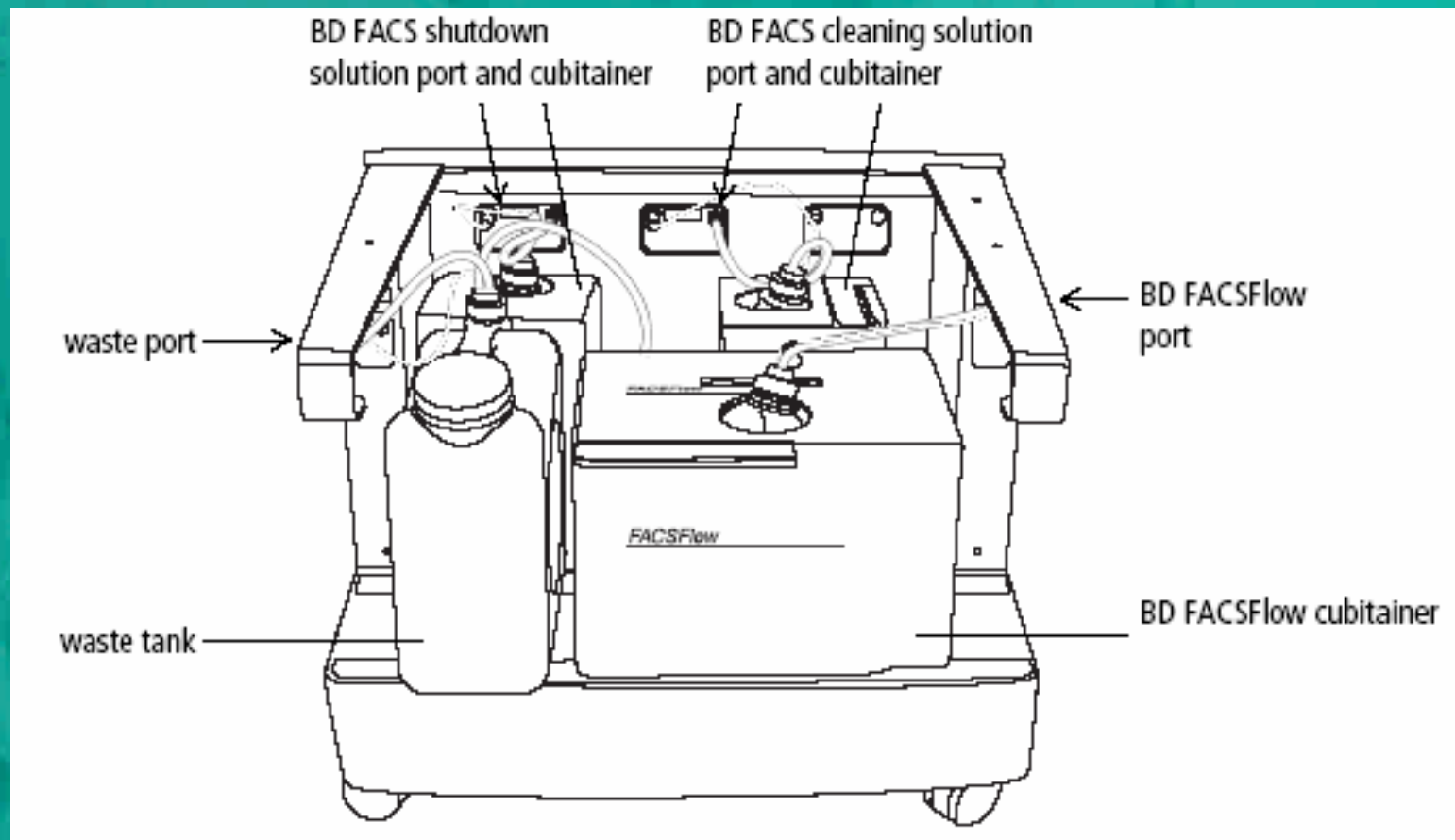
Helping all people
live healthy lives

Helping all people live healthy lives

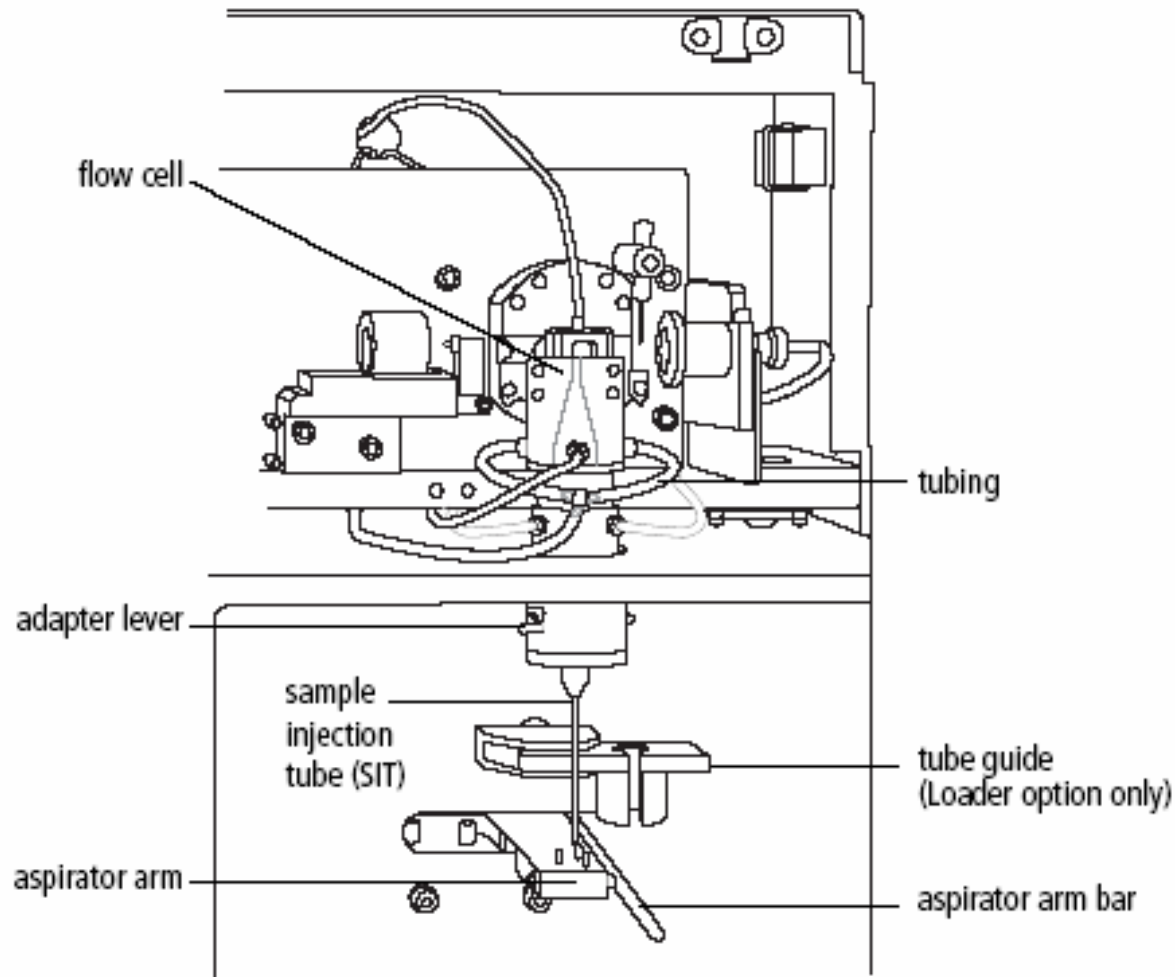
- ❑ Compressor
- ❑ 10 or 20 L sheath reservoir
- ❑ 5 L rinsing solution
- ❑ 5 L shutdown solution

Helping all people
live healthy lives

Figure 1-2 Sample injection tube

10,000 events/sec

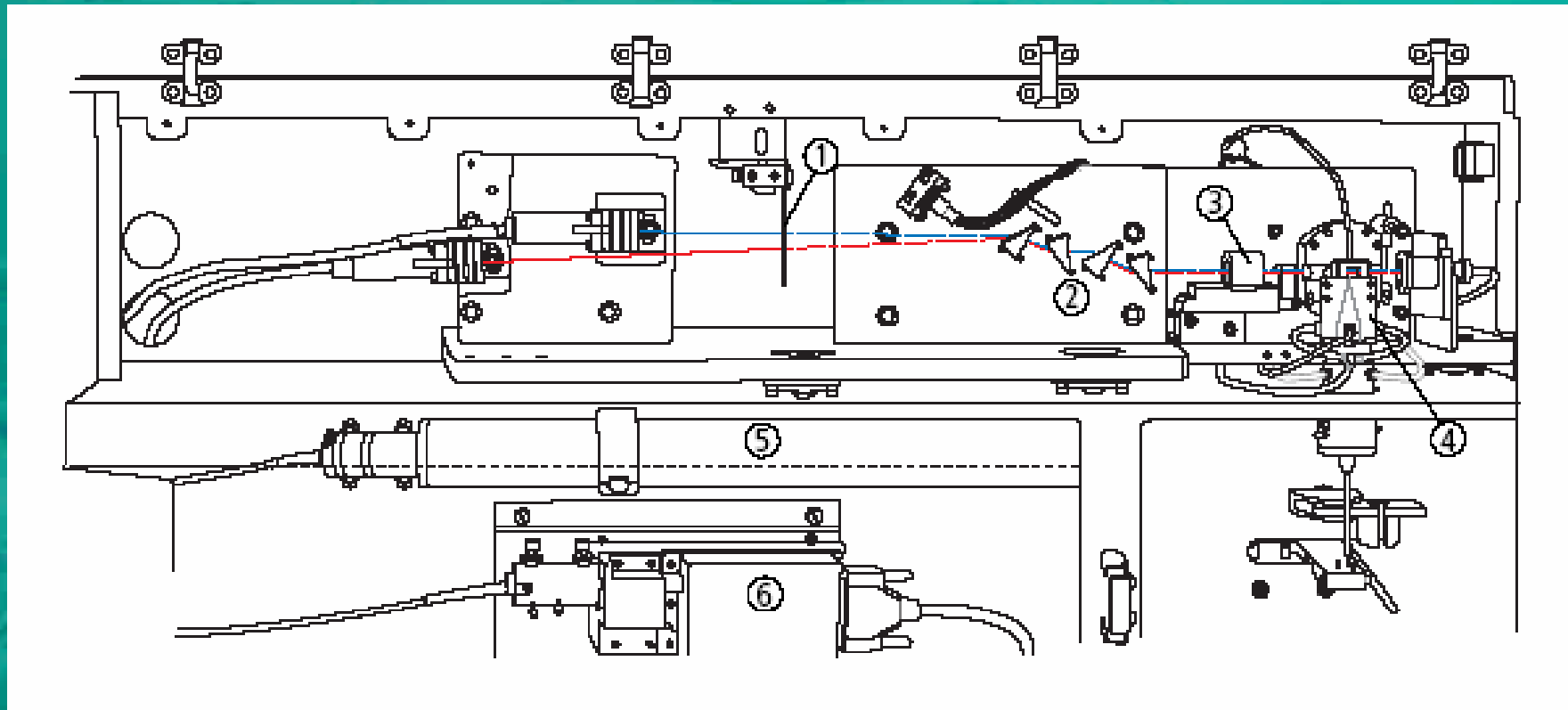
Helping all people
live healthy lives

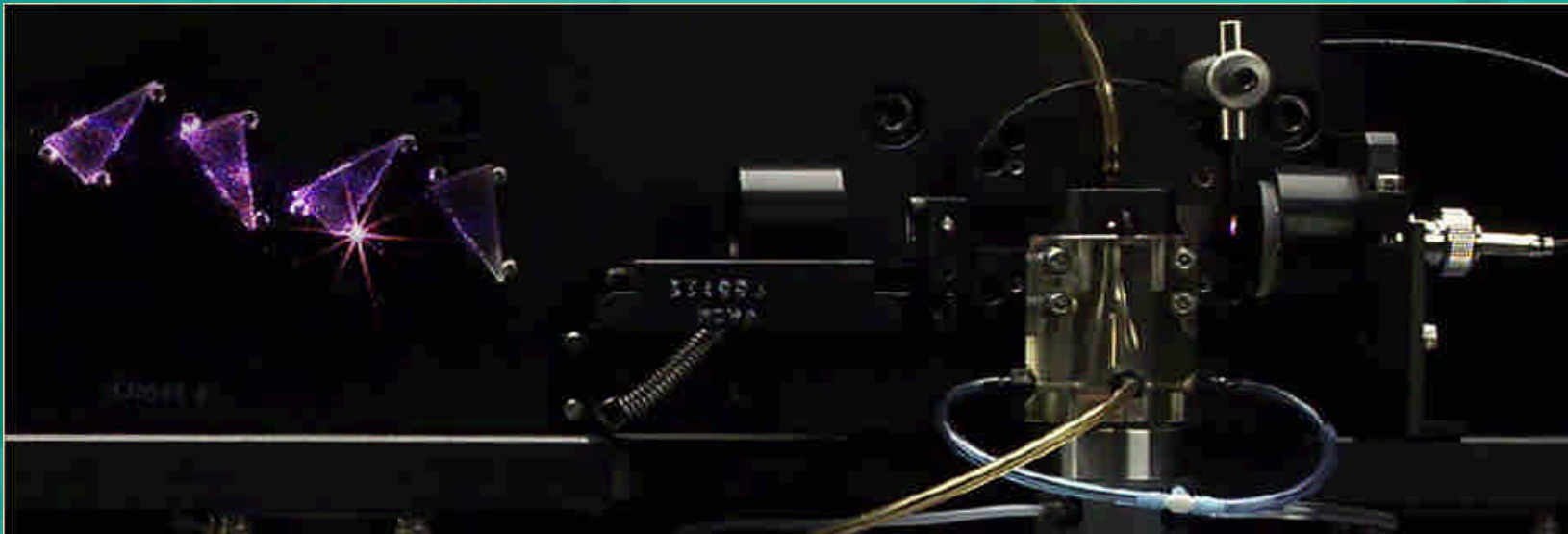
Helping all people
live healthy lives

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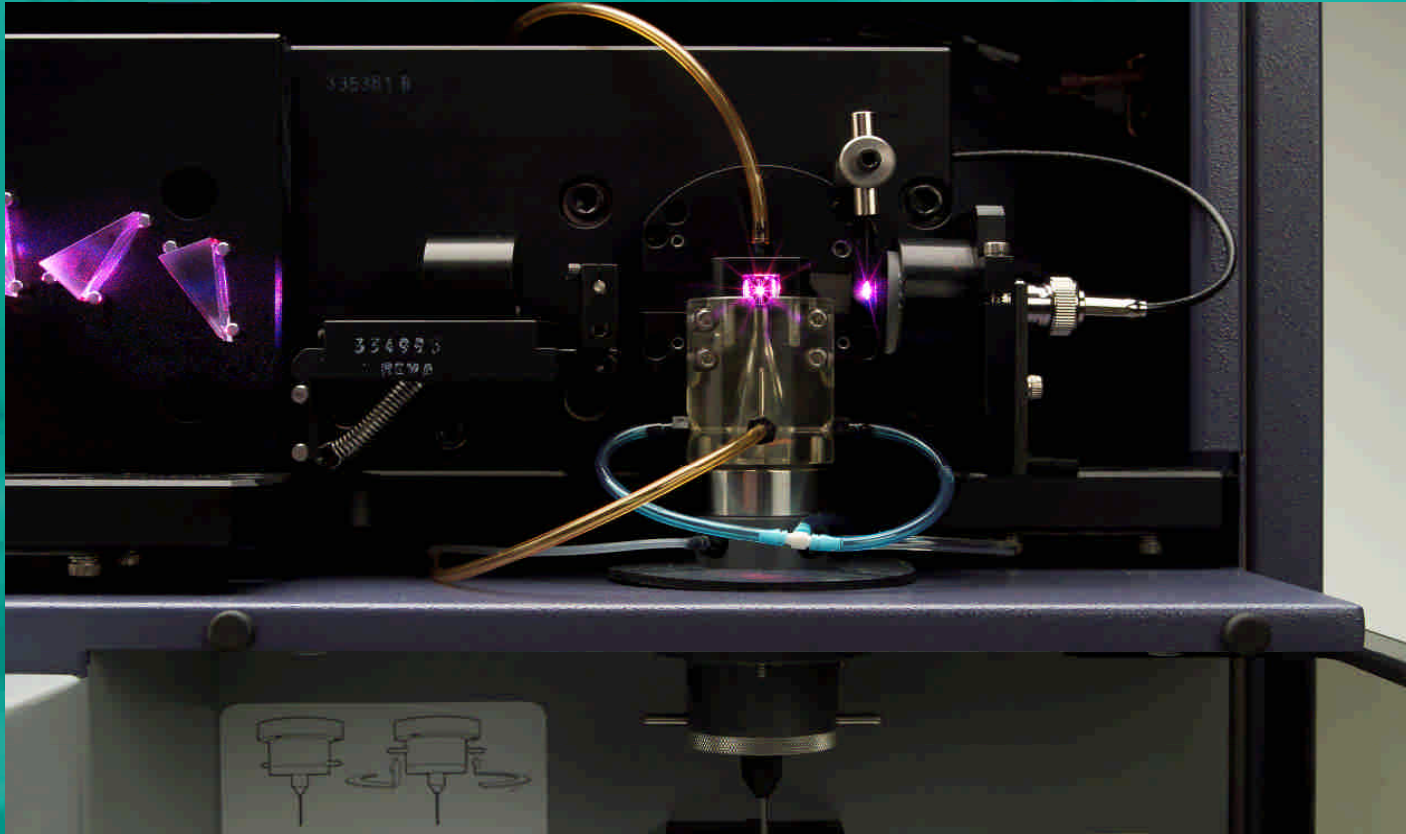
Sensitivity : 100 FITC MESF, 50 PE MESF

Helping all people
live healthy lives

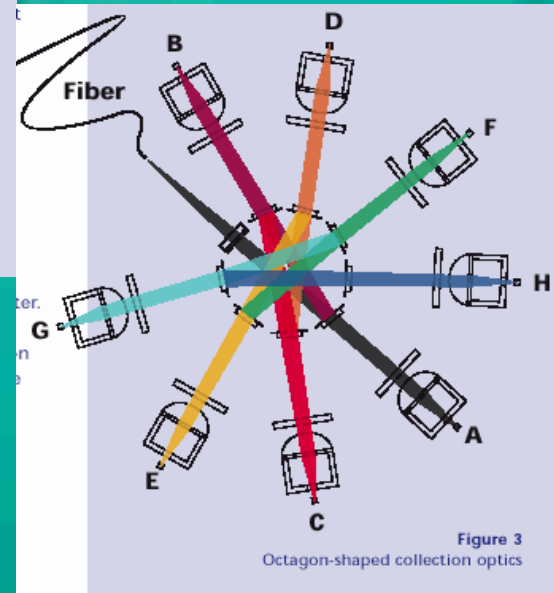
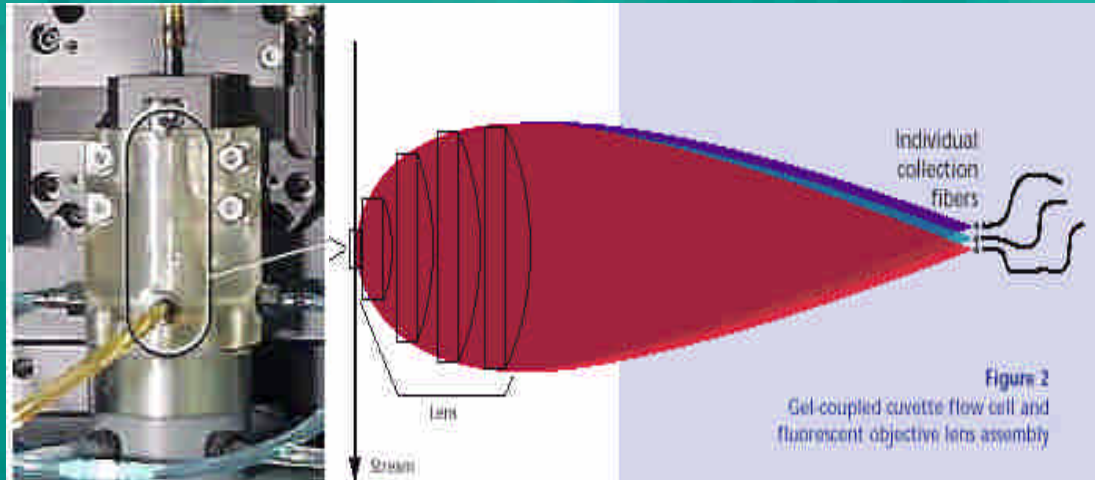
Helping all people
live healthy lives

Helping all people
live healthy lives

Cuvette Flow Cell & Fiber Optics

Helping all people
live healthy lives

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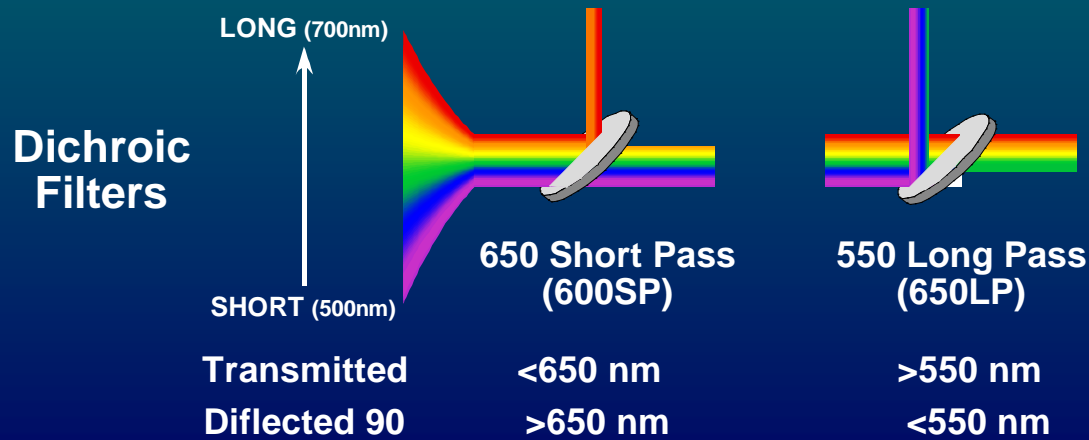
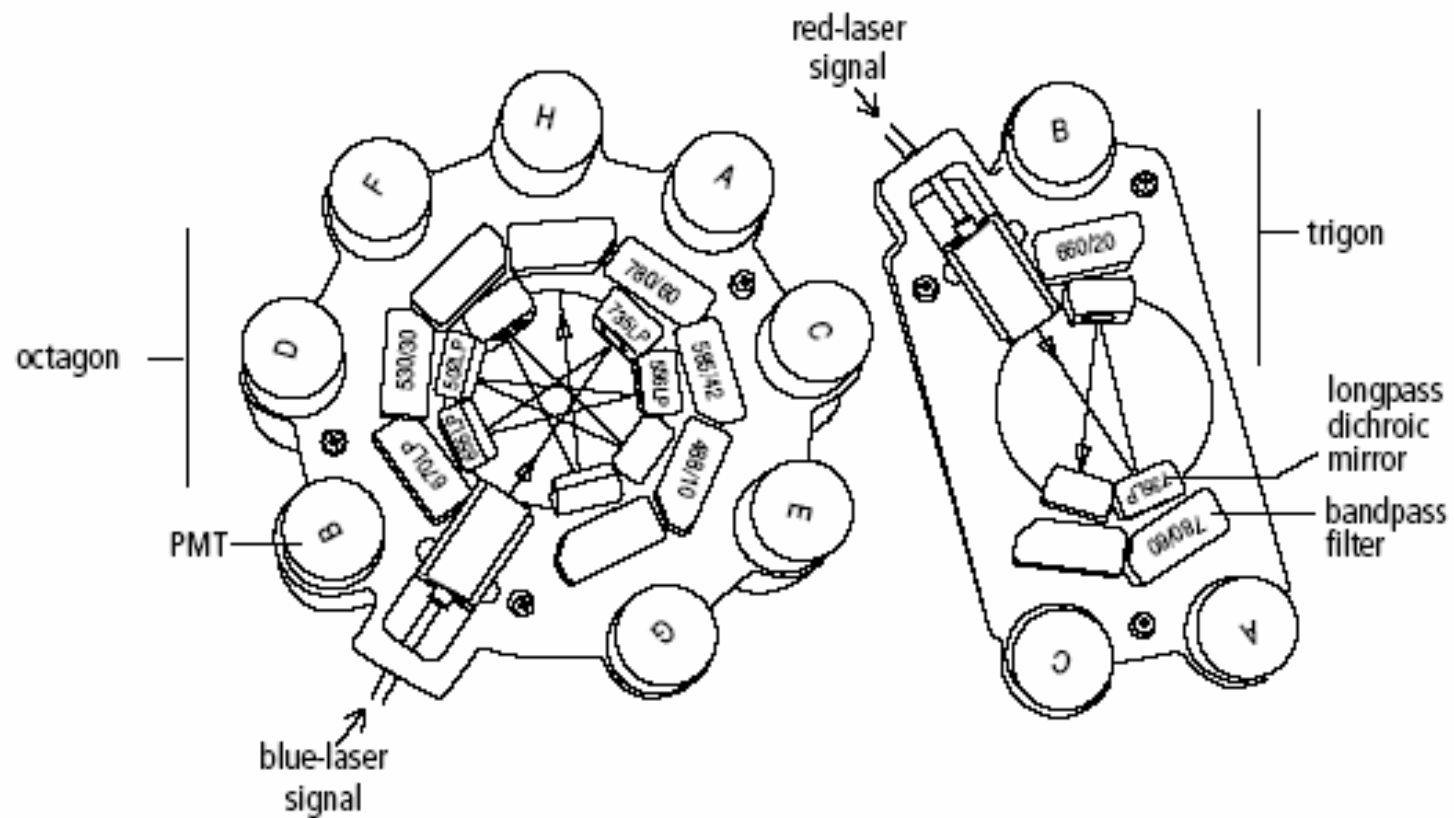
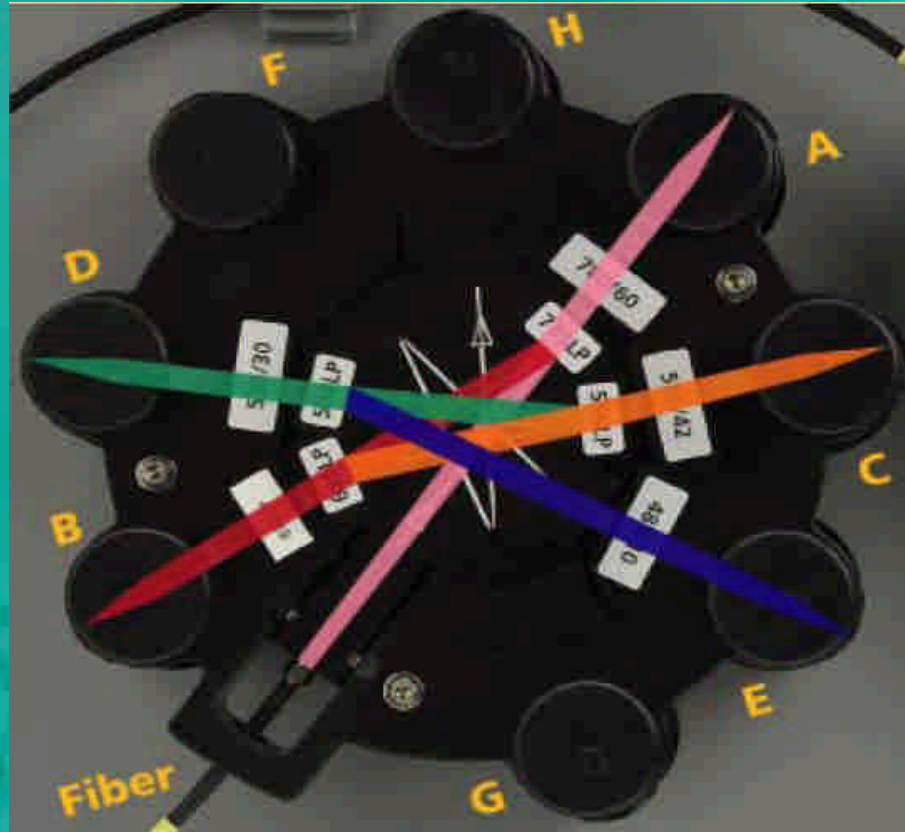
Figure 1-5 Octagon and trigon detector arrays

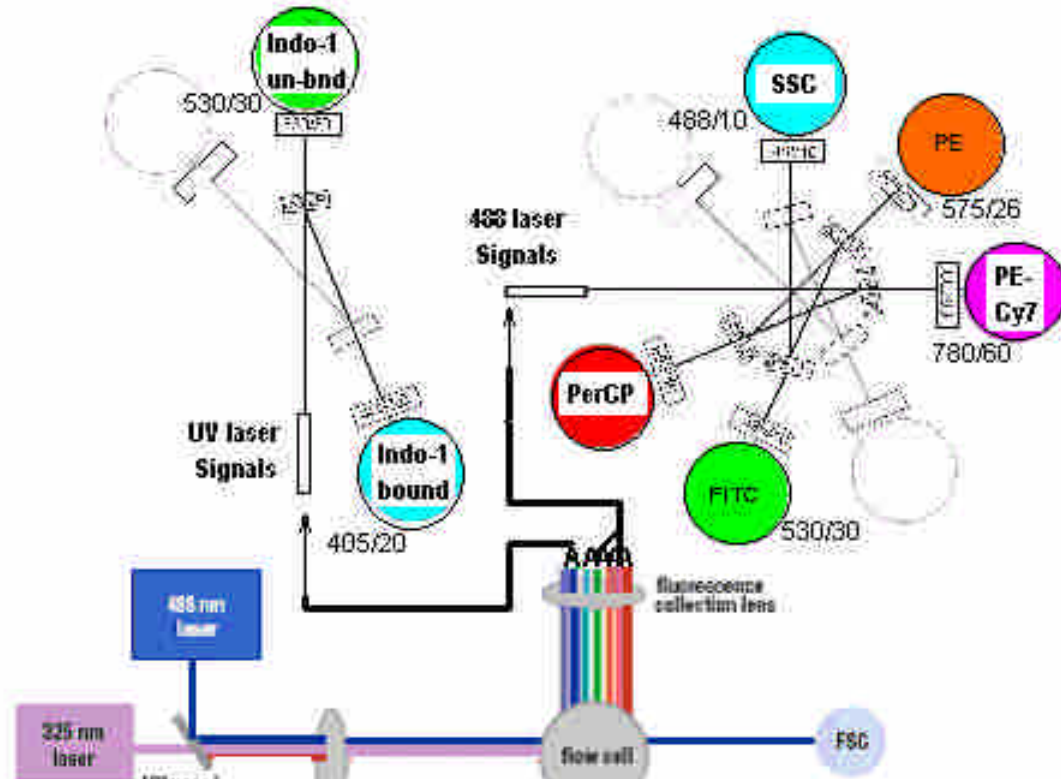
Helping all people
live healthy lives

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Helping all people
live healthy lives

Laser	Wavelength (nm)	Min. Power ^a (mW)	Commonly Used Fluorochromes
Coherent® Sapphire™ Solid State	488 (blue)	20	FITC, PE ^b , PE-Texas Red®, PerCP, PerCP-Cy5.5, PE-Cy7, PI
JDS Uniphase™ HeNe Air Cooled	633 (red)	17	APC, APC-Cy7

Helping all people
live healthy lives

BD FACSCanto – Default dyes

Instrument	Laser	Excitation Laser Line (nm)	Fluorescence Channel	Fluorochromes from BD Biosciences	
BD FACSCanto Flow Cytometer	Blue	488	Green	FITC	
			Yellow	PE	
			Red	PerCP	PerCP-Cy5.5
			Infra Red	PE-Cy7	
			Red	APC	
	Red	633	Red	APC	
			Infra Red	APC-Cy7	

Table 1

Validated fluorochrome combinations with the BD FACSCanto flow cytometer.

Helping all people live healthy lives

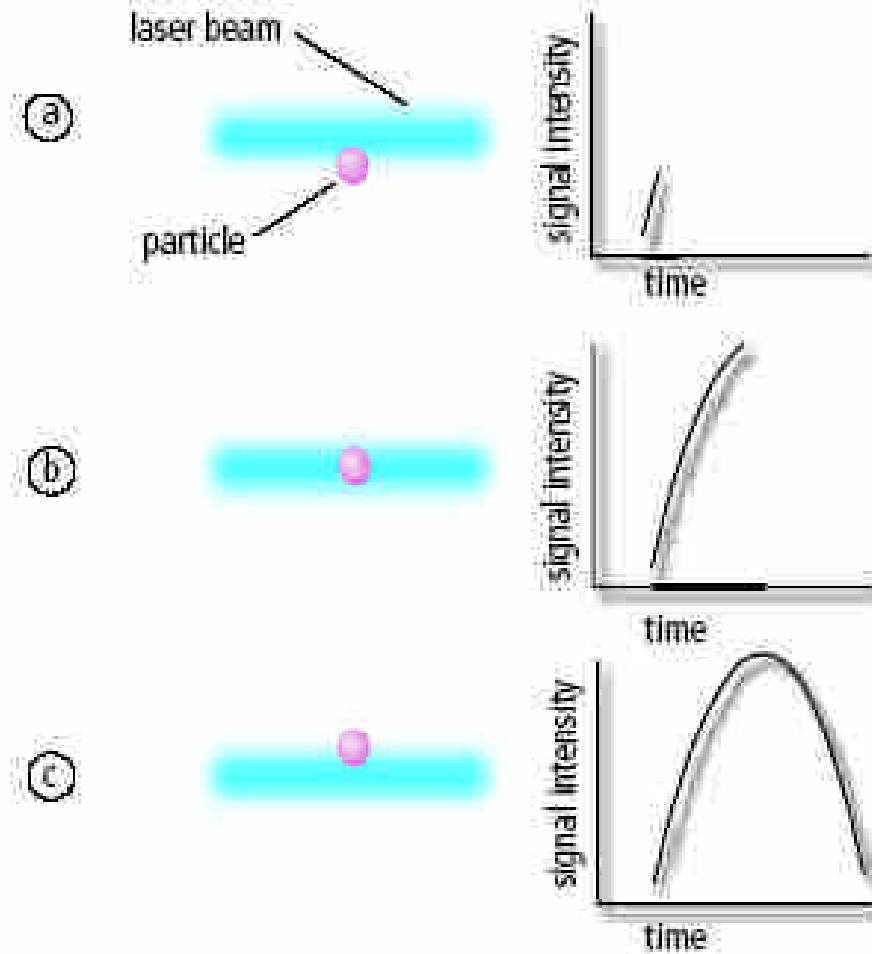
Table 1-1 Octagon and trigon optical filters

Detector Array (Laser)	PMT Position	LP Mirror	BP Filter or LP Mirror	Intended Dye
Octagon (488-nm blue laser)	A	735	780/60	PE-Cy7
	B	655	670 LP	PerCP-Cy5.5, PerCP
	C	556	585/42	PE
	D	502	530/30	FITC
	E	blank optical holder	488/10	Side scatter (SSC)
	F	blank optical holder	blank optical holder	—
	G	blank optical holder	blank optical holder	—
	H	blank optical holder	blank optical holder	—
Trigon (633-nm red laser)	A	735	780/60	APC-Cy7
	B	blank	660/20	APC

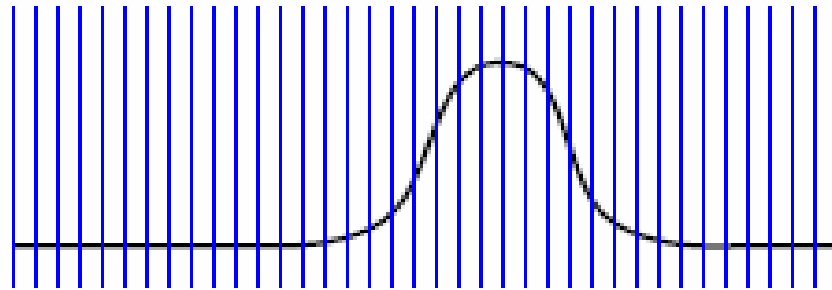
Helping all people
live healthy lives

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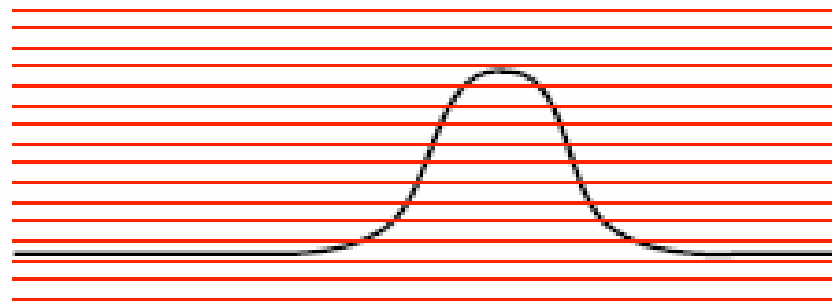
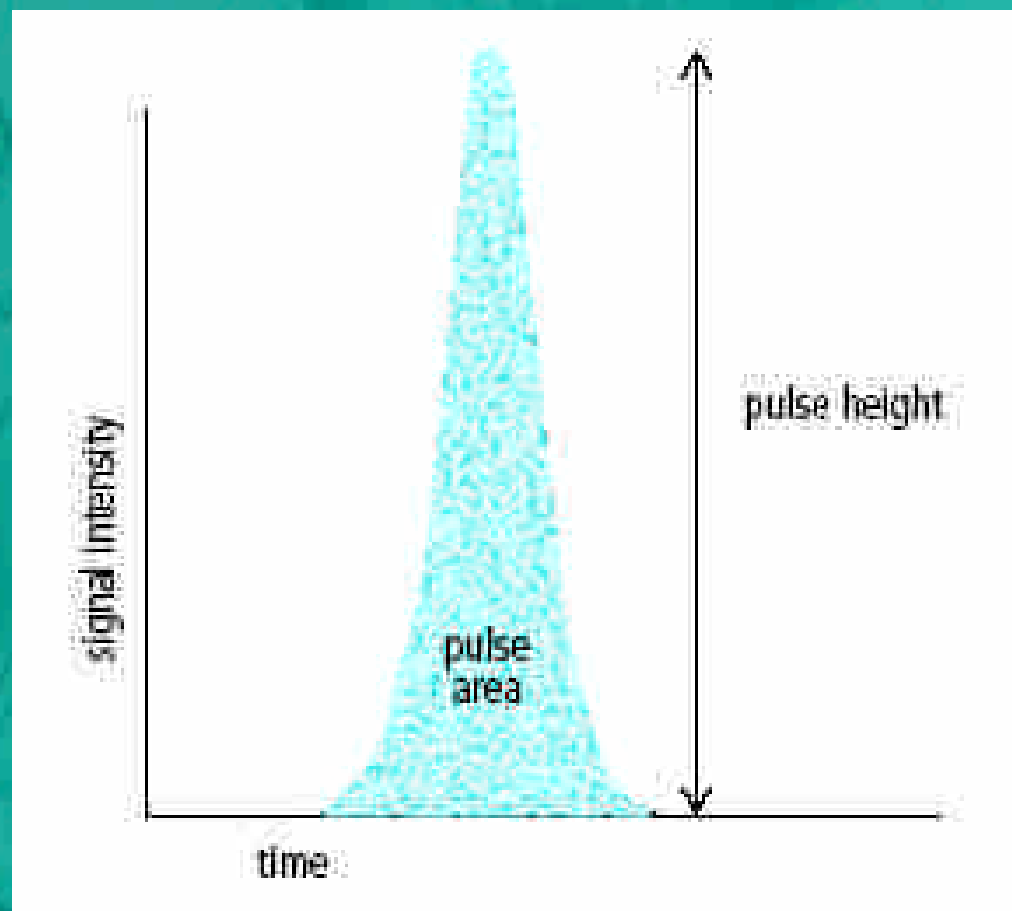
measured 10,000,000 times
every second

digitized into 16,384 levels

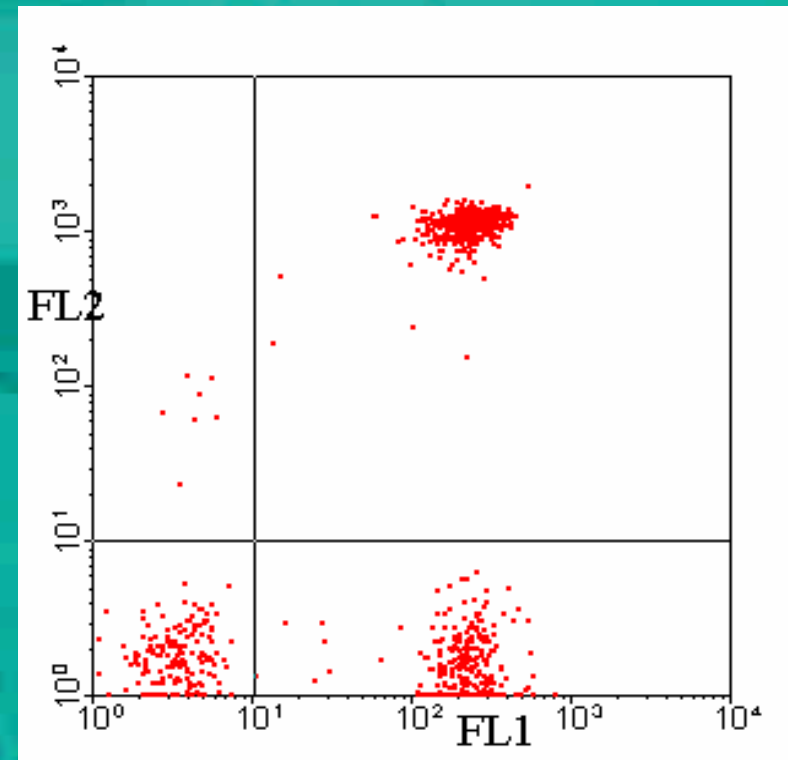
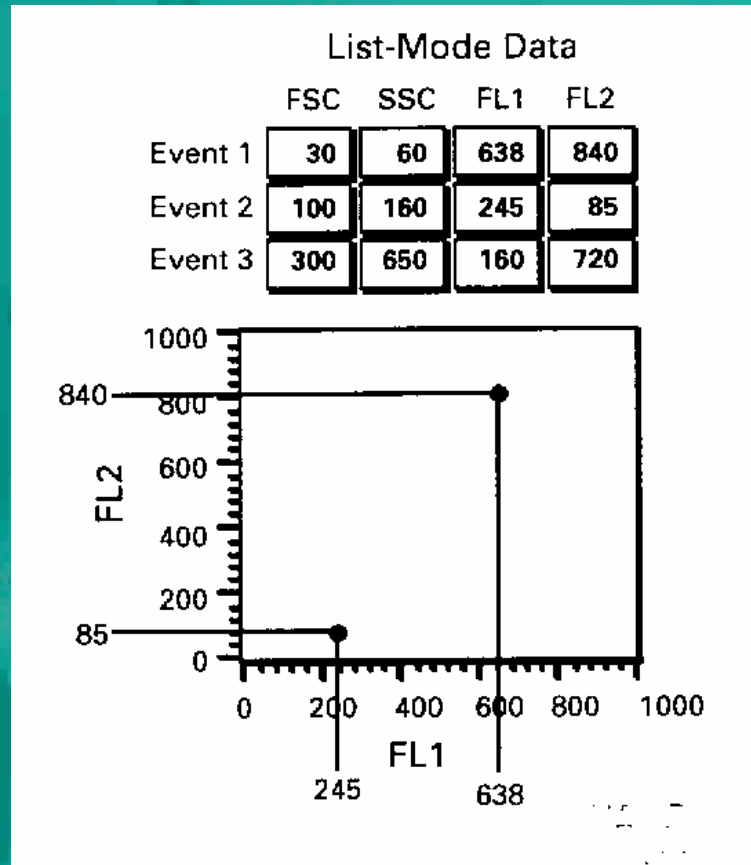
Helping all people
live healthy lives

Helping all people
live healthy lives

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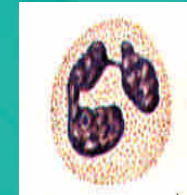
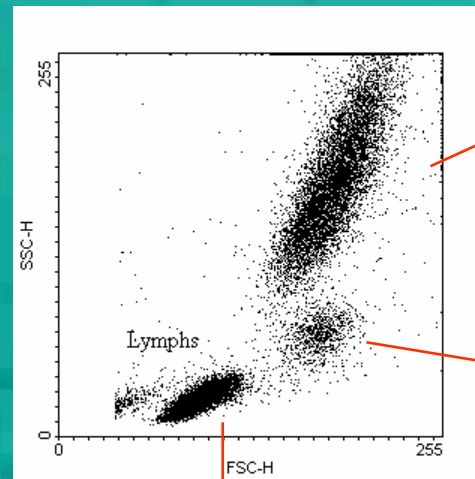
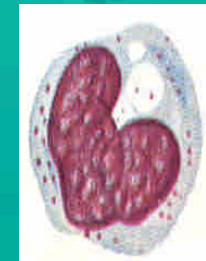
Helping all people
live healthy lives

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10 to 14 μm
Neutrophil

15 to 20 μm
Monocyte

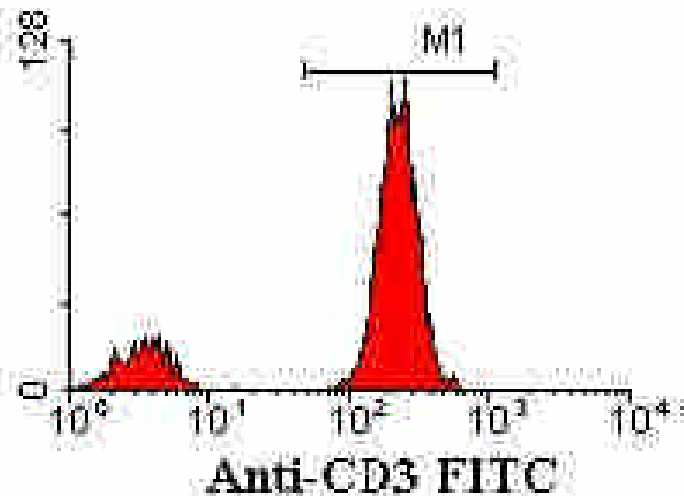
8 to 10 μm
Lymphocyte

Helping all people
live healthy lives

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File: IS20924004
Date: 24-Sep-93
Total Events 6000
Gated Events 2785

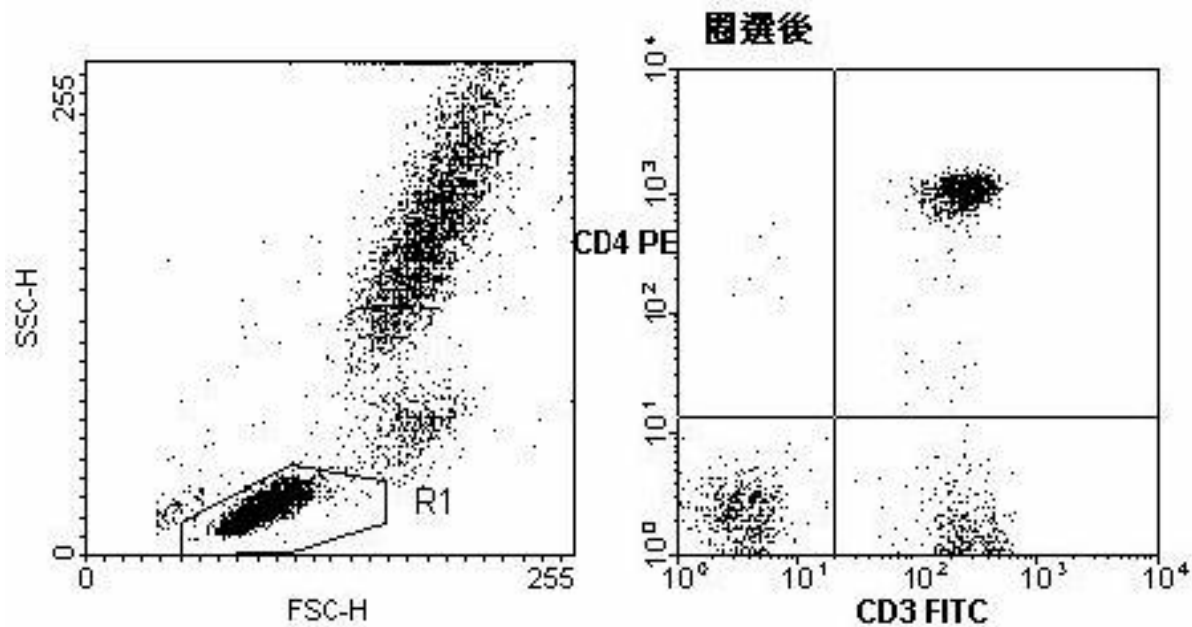
M	Low, High	Events	%Gated	GMean	CV
0	0, 255	2785	100.00	106.55	35.70
1	108, 196	2270	81.51	231.20	6.06

CV=S.D./Mean

Helping all people
live healthy lives

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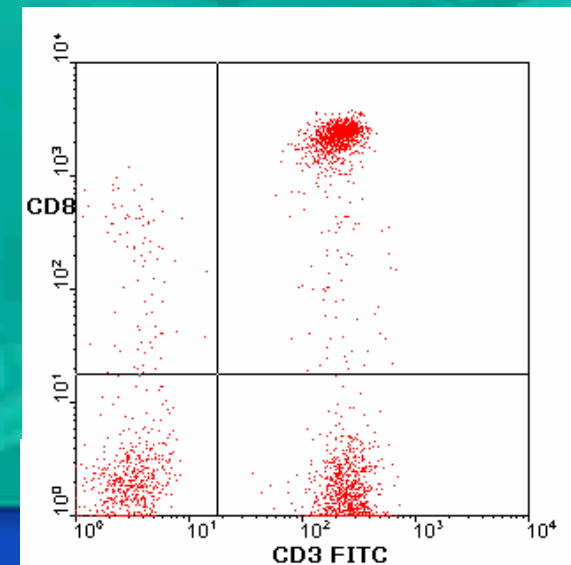
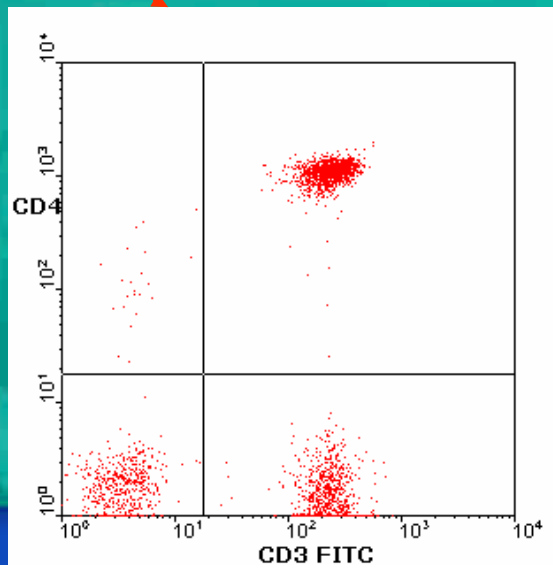
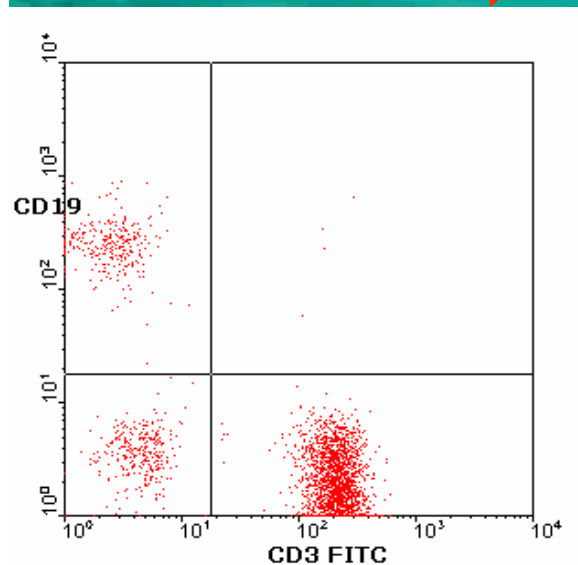
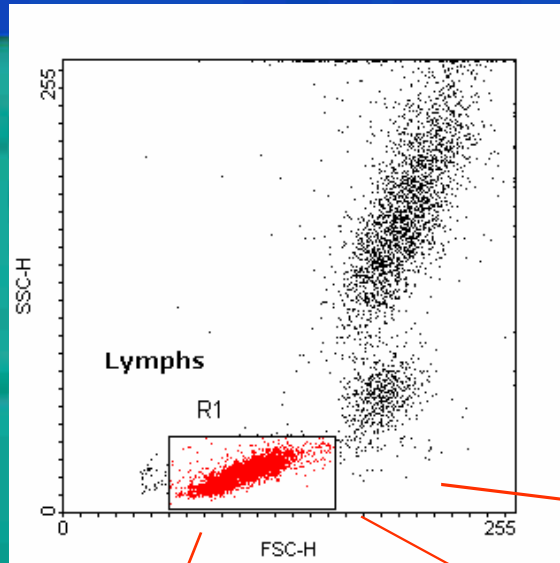
四象限統計 Quadrant Statistics

Gates: R1				
Gated Events: 2354				
Quad Stats				
FL1-H(Log) vs FL2-H(Log)				
Quadrant x,y: 83,72				
Quad	X-Mean	Y-Mean	Events	%Gated
UL	5.2	262.6	6	0.25
UR	252.9	977.9	1139	48.39
LL	3.4	2.0	508	21.58
LR	288.1	1.3	701	29.78

Helping all people
live healthy lives

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Running Samples with BD FACSDiva Software

The following topics are covered in this chapter:

- Instrument Startup on page 54
- Instrument Quality Control on page 57
- Optimization of Instrument Settings on page 68
- Data Recording and Analysis on page 79
- Daily Shutdown on page 90

Helping all people
live healthy lives

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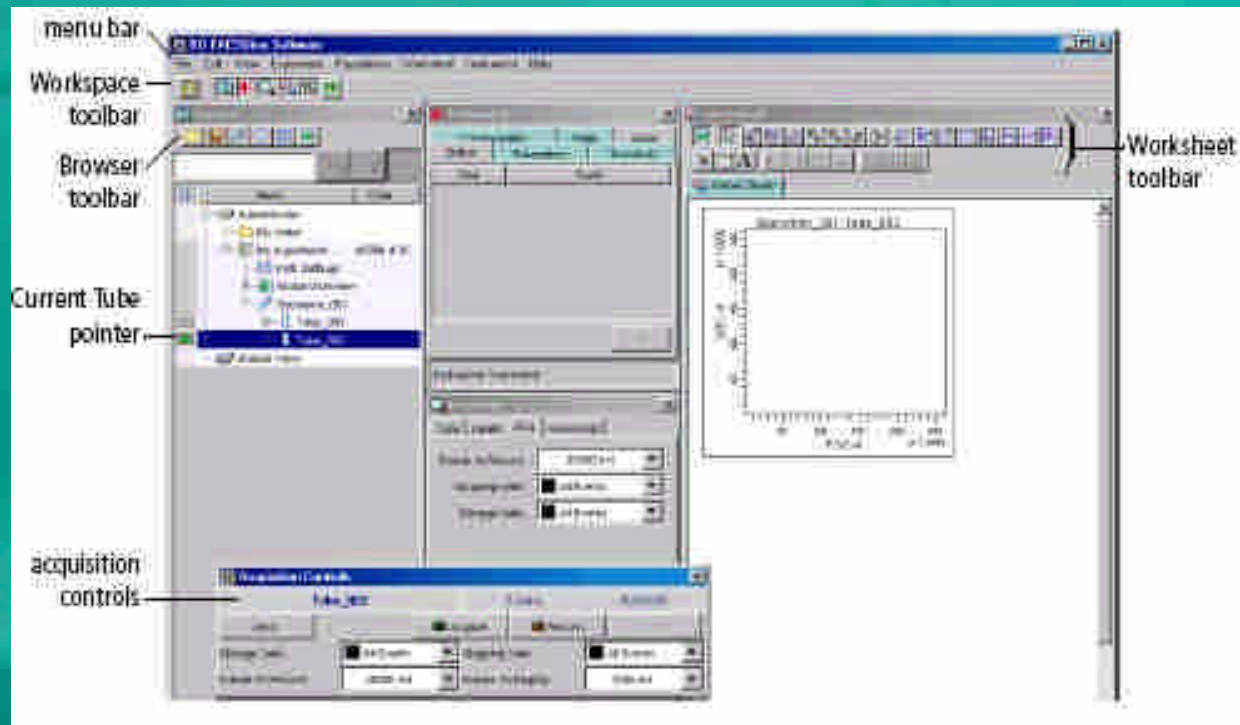
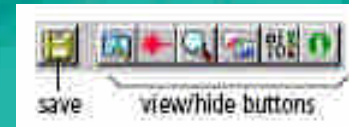
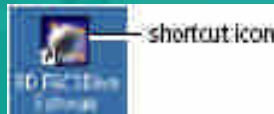
Helping all people
live healthy lives

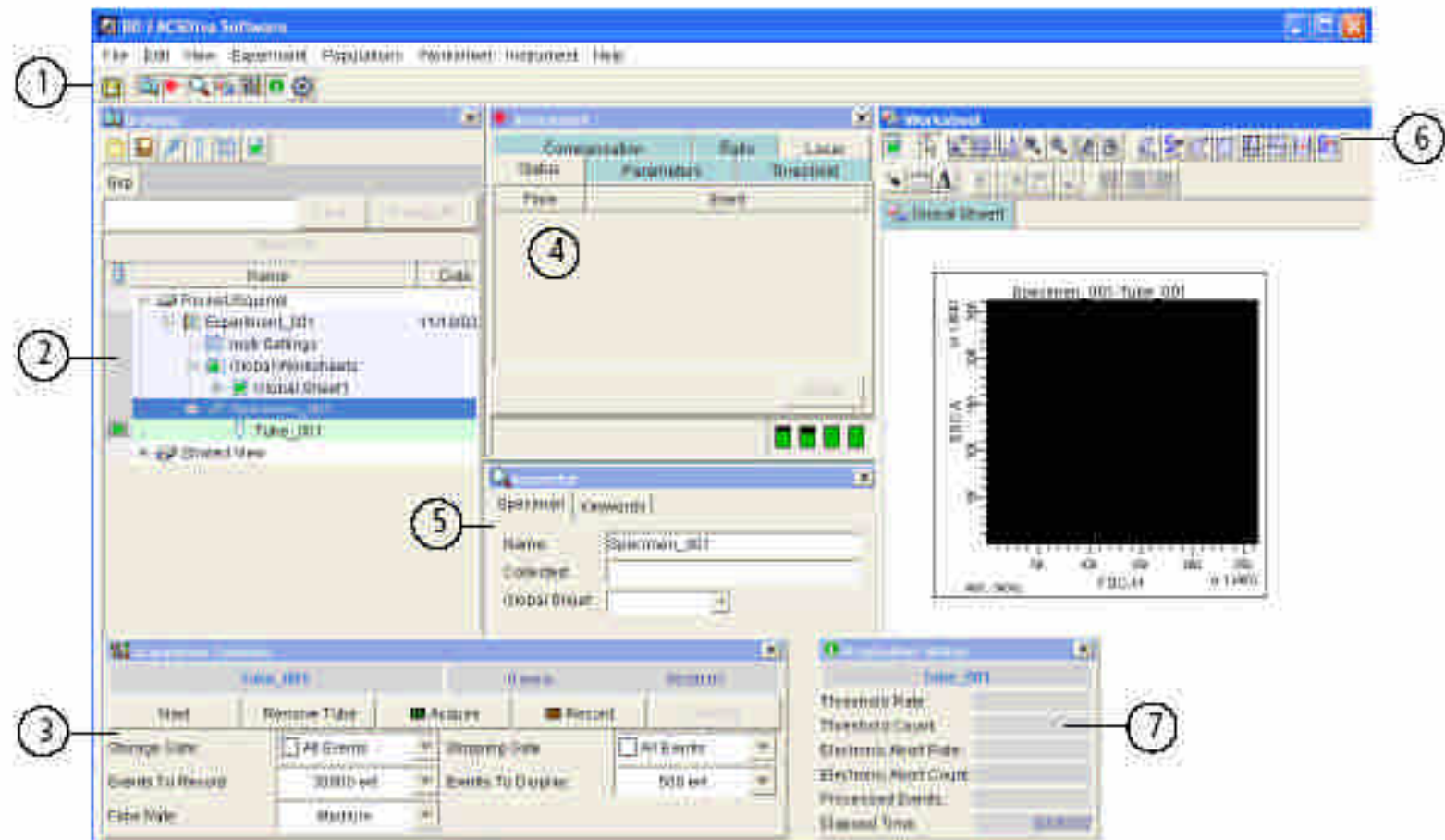
User Log In

Helping all people
live healthy lives

FACSDiva Workspace

Figure 2-1 BD FACSDiva workspace

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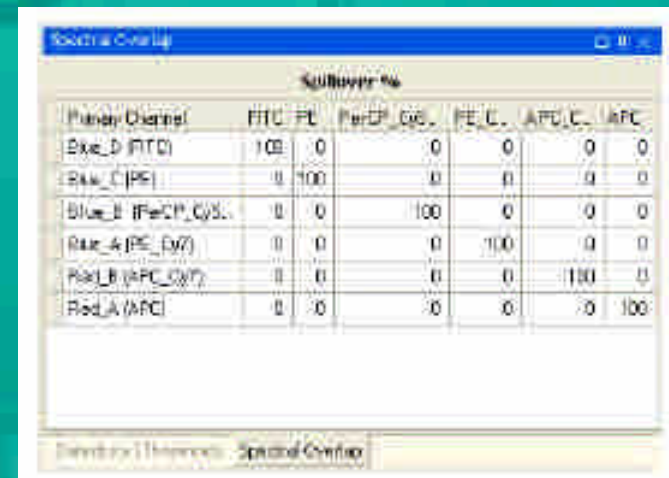
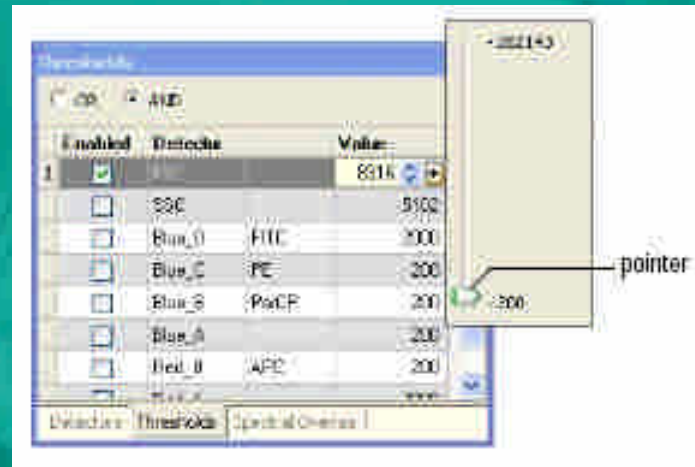
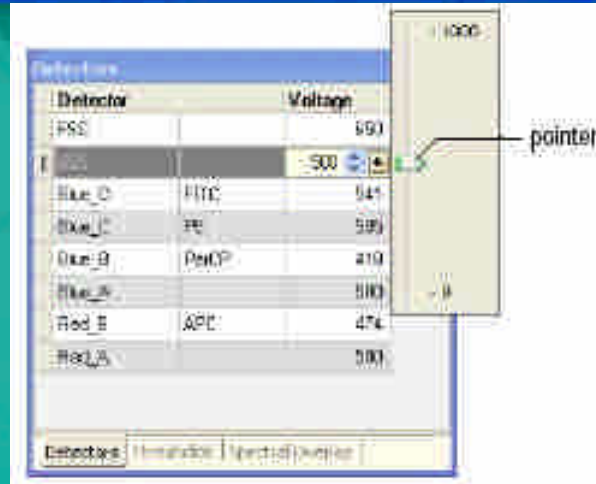
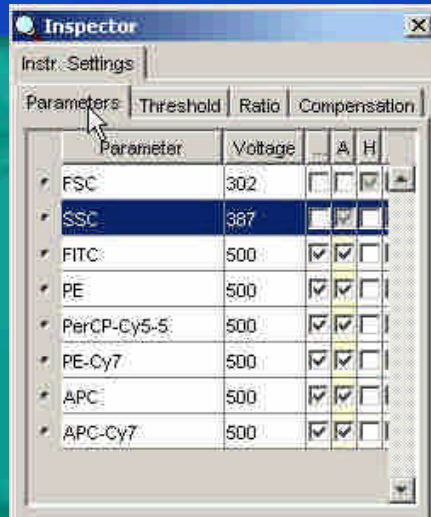
Instrument	Laser	Excitation Laser Line (nm)	Fluorescence Channel	Fluorochromes from BD Biosciences	
BD FACSCanto Flow Cytometer	Blue	488	Green	FITC	
			Yellow	PE	
			Red	PerCP	PerCP-Cy5.5
			Infra Red	PE-Cy7	
	Red	633	Red	APC	
			Infra Red	APC-Cy7	

Table 1
Validated fluorochrome combinations with the BD FACSCanto flow cytometer.

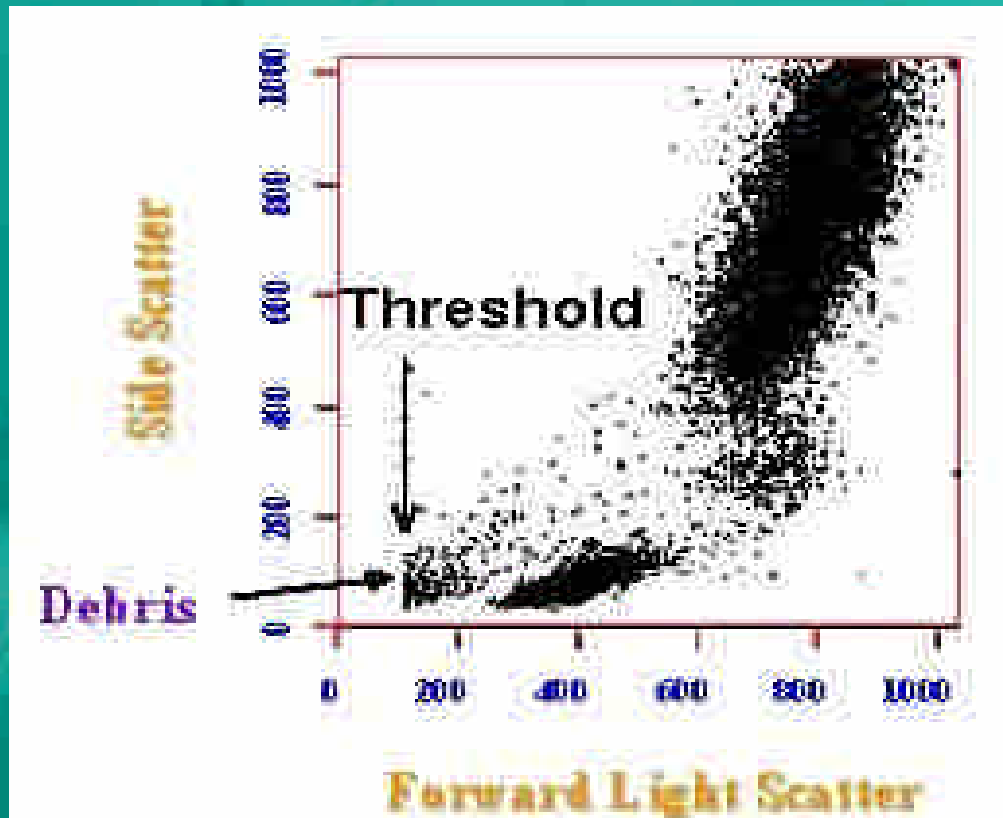
Helping all people
live healthy lives

Instrument Set Up

Helping all people live healthy lives

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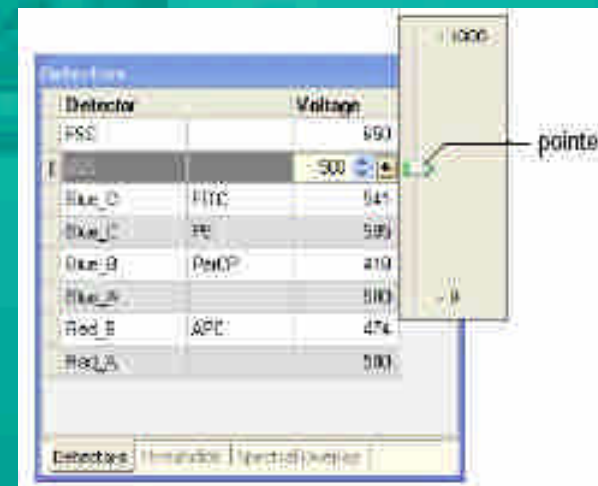
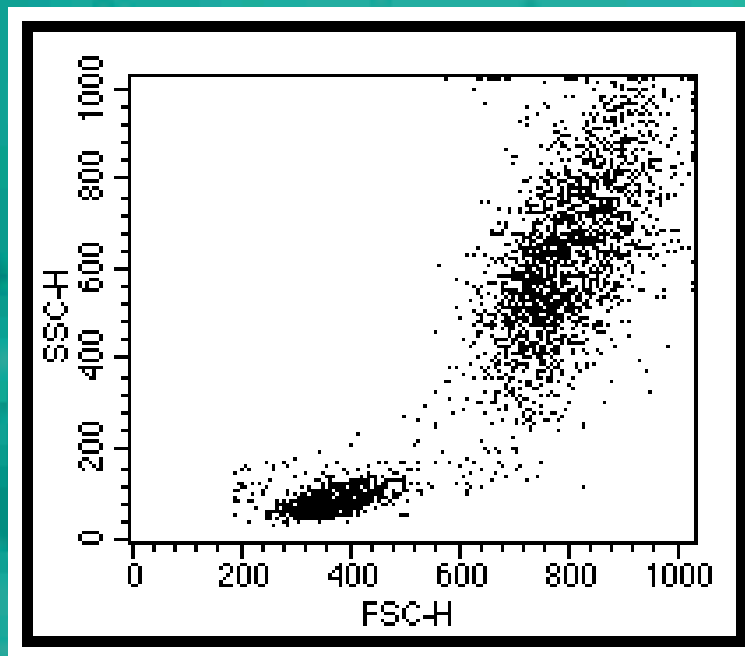
A screenshot of a software interface showing a table of parameters and a value adjustment slider. The table has columns for 'Enabled', 'Detects', and 'Value'. A slider is positioned over the 'Value' column, with a 'pointer' label pointing to the value '300'.

Enabled	Detects	Value	
<input checked="" type="checkbox"/>		831k	
<input type="checkbox"/>	SSC	5100	
<input type="checkbox"/>	Blue_1	FITC	200
<input type="checkbox"/>	Blue_C	PE	200
<input type="checkbox"/>	Blue_B	PerCP	200
<input type="checkbox"/>	Blue_A		200
<input type="checkbox"/>	Red_1	APC	200
<input type="checkbox"/>	Red_2		200

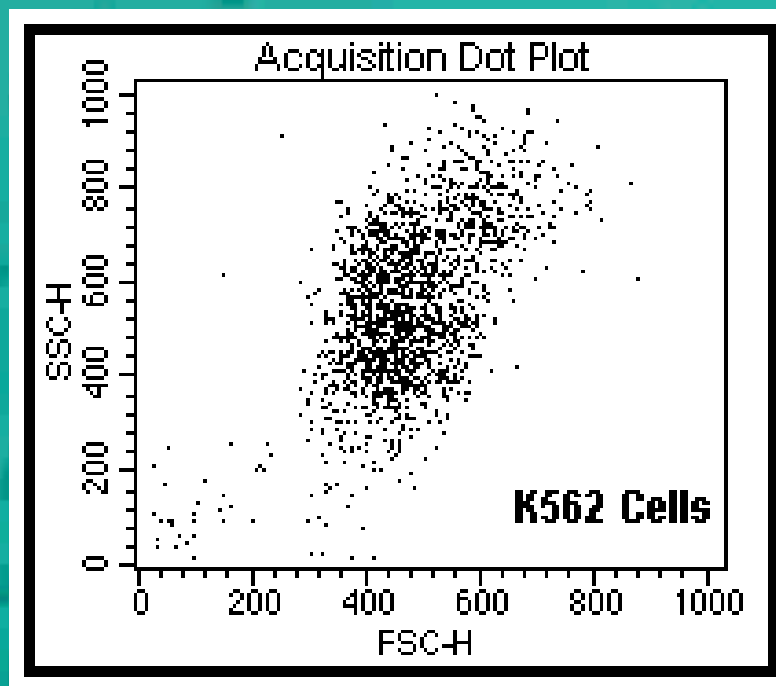
Helping all people live healthy lives

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Helping all people
live healthy lives

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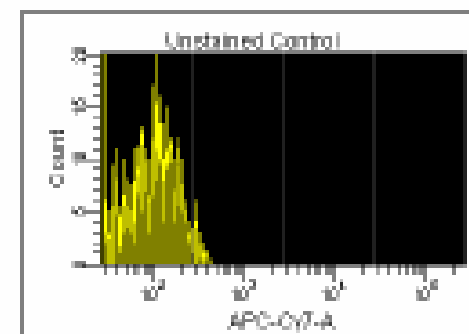
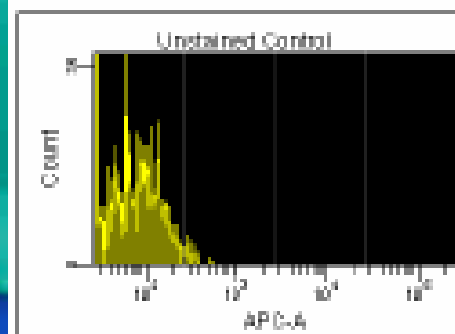
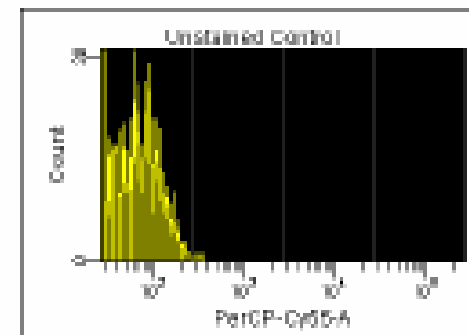
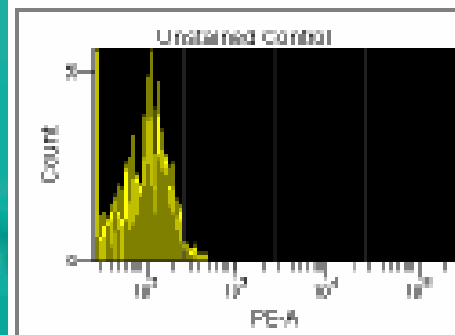
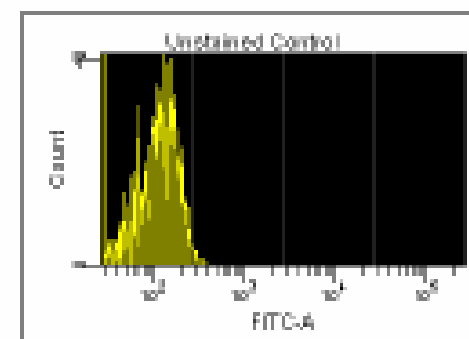
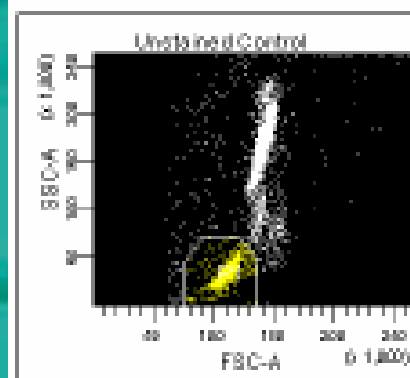
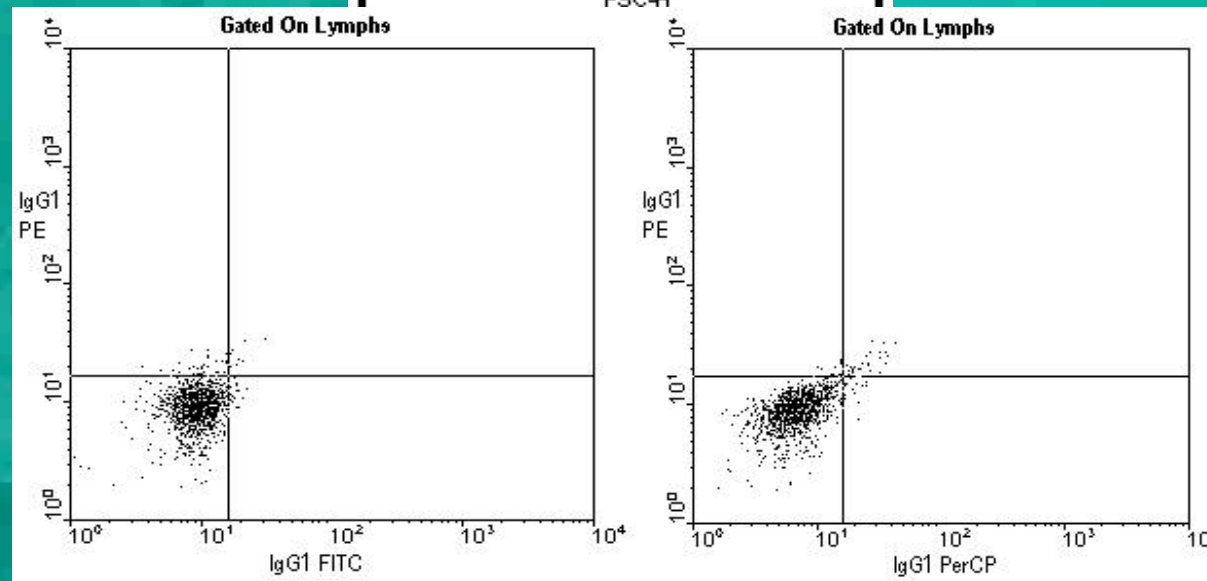
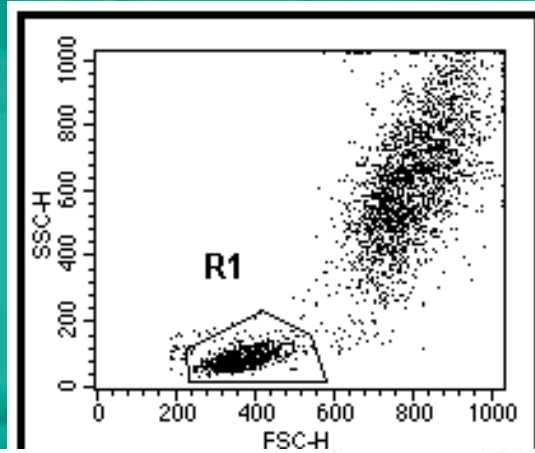
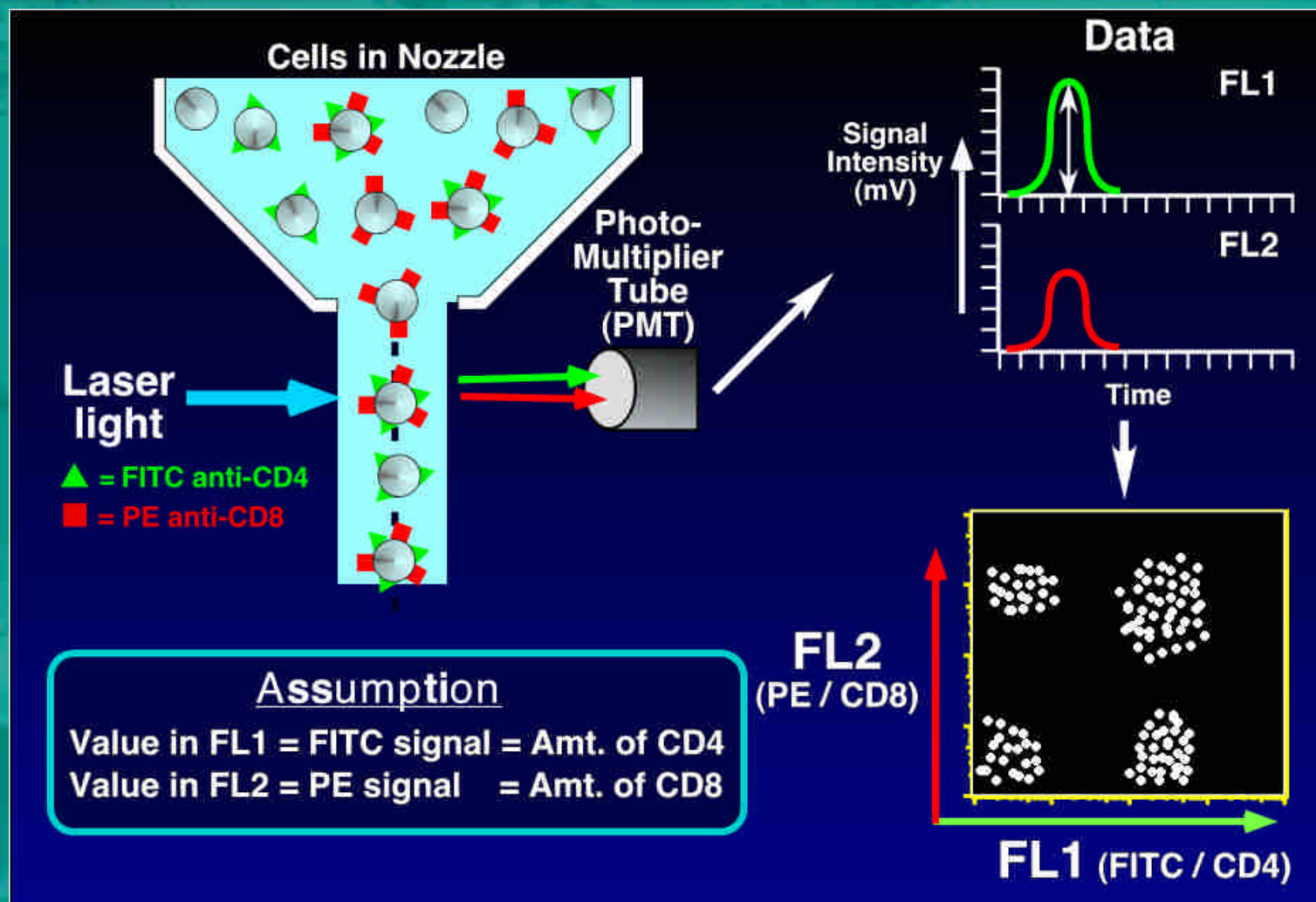
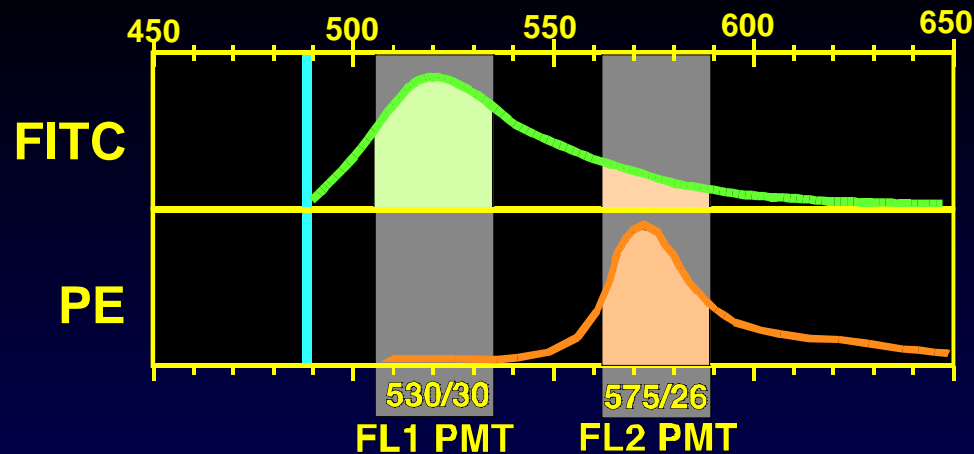
$$FL1 = FITC + x\% PE$$

$$FL2 = PE + y\% FITC$$

Remember the basic assumption of flow analysis:

The signal in FL1= the signal from FITC and only FITC and
the signal in FL2= the signal from PE and only PE.

This is NOT TRUE for the raw data!

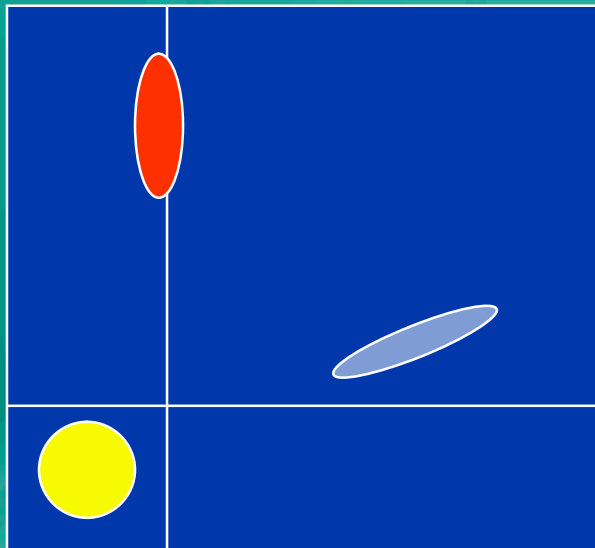
The process by which each fluorescence channel is "corrected" for this spectral overlap is termed *Fluorescence Compensation*

Helping all people
live healthy lives

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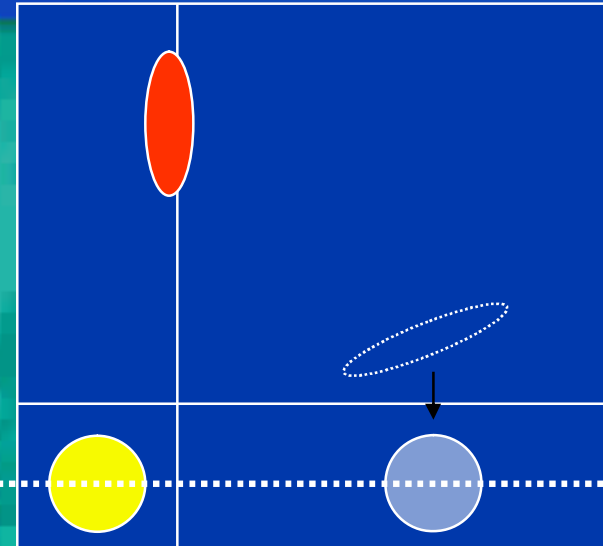
FL-2(PE)

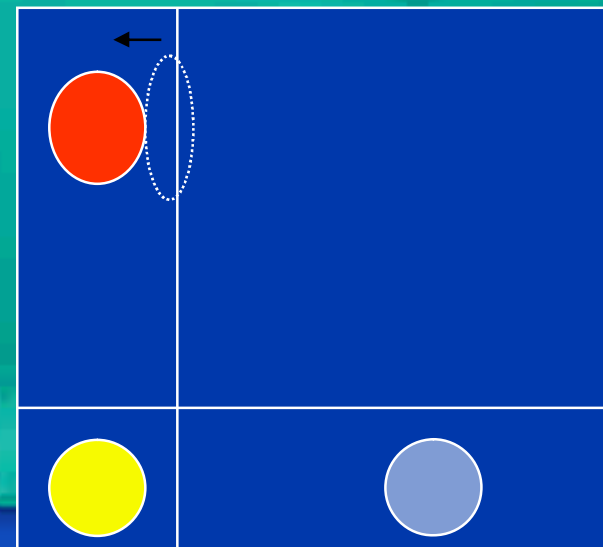
FL-1(FITC)

FL2 - %FL1
(18 - 30%)

FL1 - %FL2
(0 - 1%)

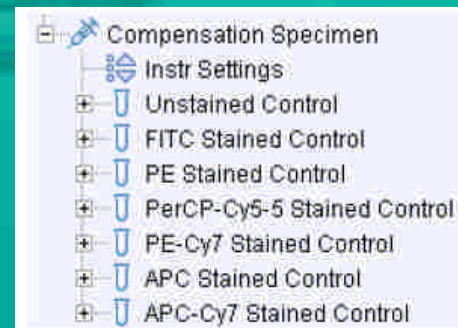
Create Compensation Controls

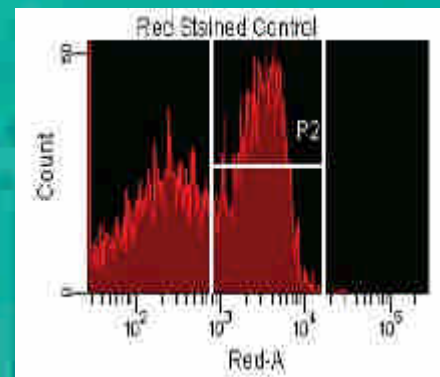
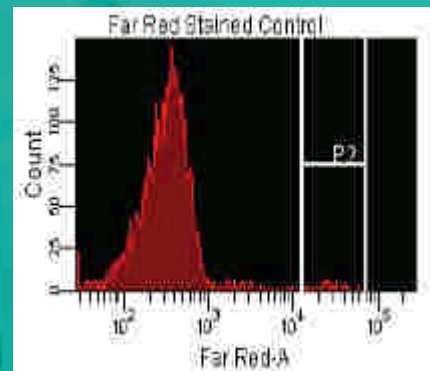
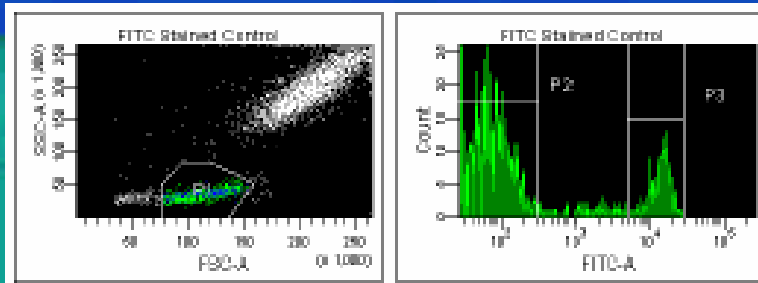
Include separate unstained control tube/well

Fluorophore	Label
FITC	Generic
PE	Generic
PerCP-Cy5-5	Generic
PE-Cy7	Generic
APC	Generic
APC-Cy7	Generic

Add Delete OK Cancel

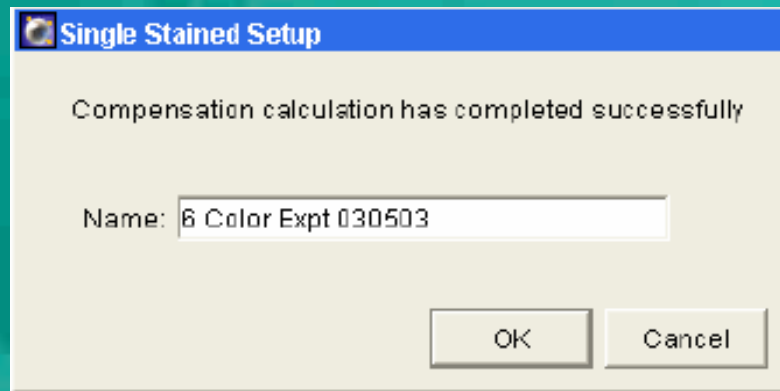
Helping all people
live healthy lives

Helping all people
live healthy lives

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A window titled "Spectral Overlay" with a blue header bar. It displays a table with the following data:

Primary Channel	FTC	FL	PerCP_Cy5	PE_C	APC_C	APC
Exc_D (FTC)	100	0	0	0	0	0
Exc_C (PE)	0	100	0	0	0	0
Blue_C (PerCP_Cy5)	0	0	100	0	0	0
Red_A (PE_Cy7)	0	0	0	100	0	0
Red_B (APC_Cy7)	0	0	0	0	100	0
Red_A (APC)	0	0	0	0	0	100

At the bottom of the window, there is a dropdown menu currently set to "Spectral Overlay".

Helping all people
live healthy lives

FACSCanto 各部構造

Helping all people
live healthy lives

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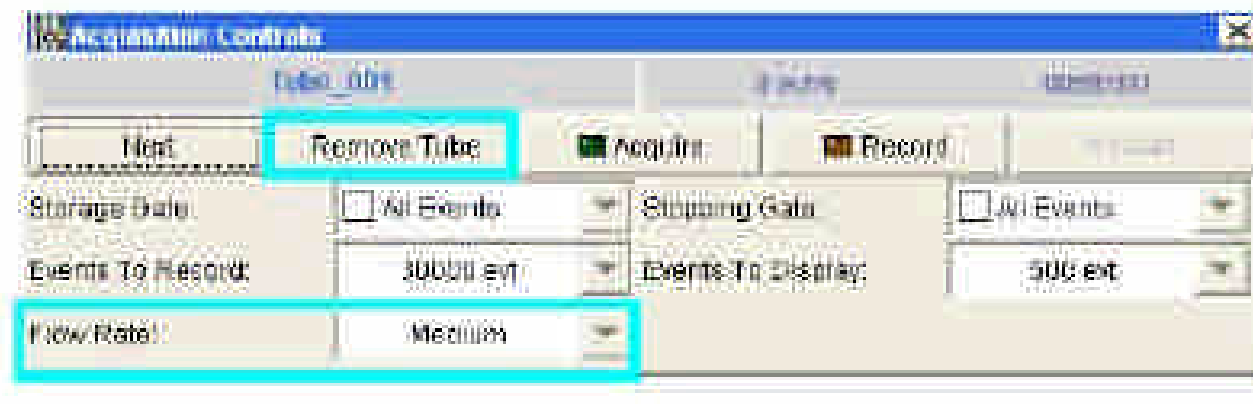
- Fisher Hematology Diluent
- Isoton III
- Isolac D
- DI water

Helping all people
live healthy lives

Figure 2-7: Acquisition controls unique to the BD FACSCanto

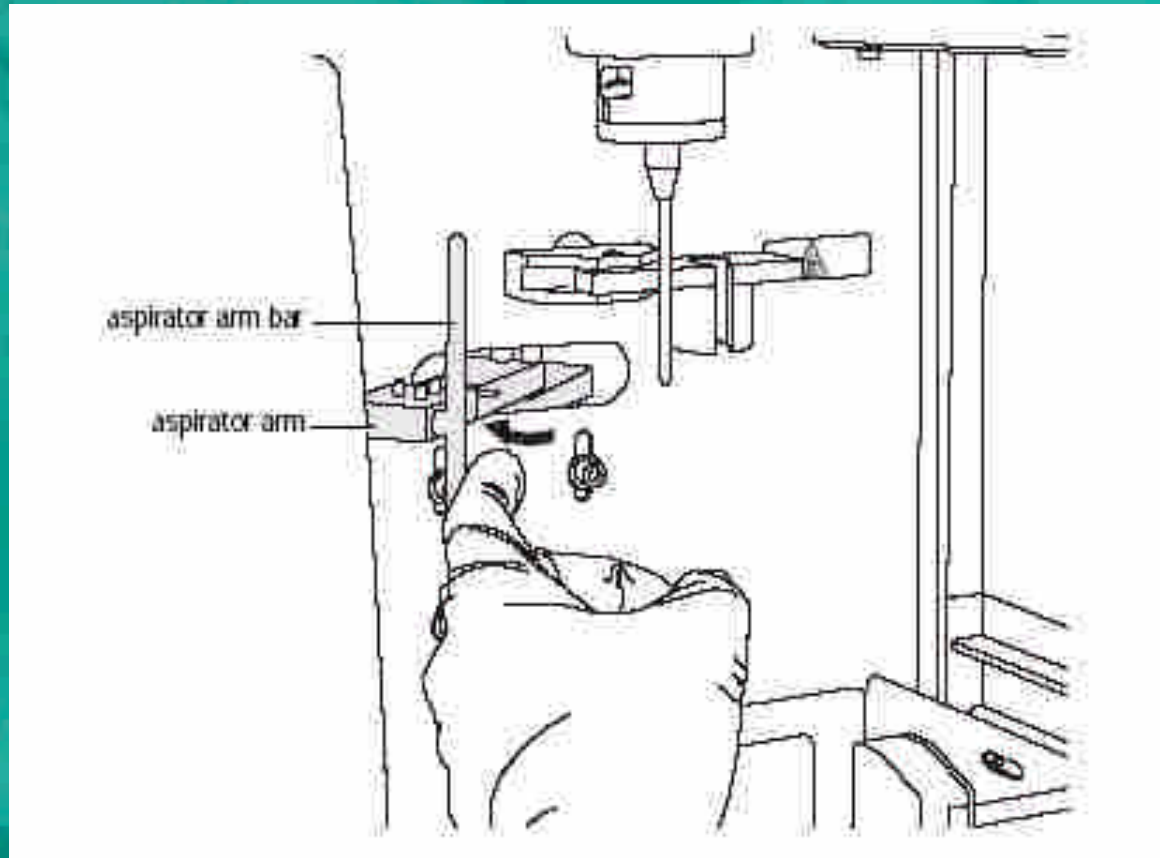
Low: 样品流速= 10 μl /min
Medium: 样品流速= 60 μl /min
High: 样品流速=120 μl /min

Helping all people
live healthy lives

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Helping all people
live healthy lives

Remove Tube 正確步驟

步驟要確實：〔 to minimize carry over between samples 〕

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